

# VLP Factory™

## Virus Like Particle Expression in Insect Cells



### User Manual

*Version 1.2*

*December 2019*

---

We thank Dr. Imre Berger and his team from Bristol, for their input and assistance in preparing this manual.

## TABLE OF CONTENTS

<b>A. VLP Factory™ Kit Contents</b>	<b>- 2 -</b>
Reagents to be supplied by user	- 3 -
<b>B. Overview</b>	<b>- 4 -</b>
<b>C. New Baculovirus Tools for Virus Like Particle (VLP) Applications</b>	<b>- 10 -</b>
C.1. Transfer vectors: the Acceptor-Donor recombineering system.	- 10 -
C.2. Generating multigene expression cassettes	- 13 -
C.2.1. Multigene construction via homing endonuclease/BstXI multiplication	- 13 -
C.2.2. Multigene construction using Cre-Lox recombination	- 16 -
C.2.3. Combining HE/BstXI cycling and Cre-Lox recombination	- 17 -
C.3. Baculovirus engineered for improved protein production.	- 18 -
C.4. Introducing additional control elements	- 20 -
<b>D. Protocols</b>	<b>- 21 -</b>
D.0 Introductory remarks	- 21 -
D.1 Cloning into pACEBac or pIDx transfer vectors	- 21 -
D.2 Multiplication by using the HE and BstXI sites	- 21 -
D2.1 Protocol 1. Multiplication by using homing endonuclease/BstXI.	- 22 -
D.3 Cre-LoxP reaction of Acceptors and Donors	- 25 -
D 3.1. Protocol 2: Cre-LoxP fusion of Acceptors and Donors	- 25 -
D 3.2. Protocol 3. Deconstruction of fusion vectors by Cre	- 28 -
D.4. Transposition protocol for pACEBac derivatives (electroporation).	- 31 -
D.5. Bacmid preparation and infection of insect cells.	- 31 -
D.6. Determining baculoviral expression	- 31 -
<b>D.7. Protocol for VLP purification</b>	<b>- 31 -</b>
<b>E. Appendix</b>	<b>- 34 -</b>
E.1. Verifying bacmid integrity in DH10VLPFactory™ cells	- 34 -
E.2. Preparing competent DH10VLPFactory™ cells	- 34 -

E.2.1 Protocol for electrocompetent cells	- 34 -
E.2.2 Protocol for chemically competent cells	- 35 -
E.3. Preparing bacterial stocks from agar stabs	- 36 -
E.3. VLP Factory™ vectors: maps, sequences, restriction	- 38 -
<b>E.3.1 Acceptor vectors</b>	- 39 -
E.3.1.1 pACEBac1: 2904 bp	- 39 -
E.3.1.1 pACEBac2: 2761 bp	- 41 -
<b>E.3.2 Donor vectors</b>	- 43 -
E.3.2.1 pIDC: 2312 bp	- 43 -
E.3.2.2 pIDK: 2281 bp	- 45 -
E.3.2.3 pIDS: 2231 bp	- 47 -
E.3.3. Analytical restriction digest patterns	- 49 -
E.3.4 VLP Factory™ baculoviral genome	- 50 -
<b>F. References</b>	- 53 -
<b>G. Purchaser Notification</b>	- 55 -

## VLP Factory™ Kit Contents:

- **Plasmid acceptor vectors**

pACEBac1, pACEBac2; approx. 5 µg DNA per vial (in buffer solution)

keep at 4°C for short-term storage and in a freezer at -20°C or lower for medium- and long-term storage (take care to avoid repeated freeze-thaw cycles, e.g. by aliquotting DNA prior to freezing)

- **Plasmid donor vectors**

pIDC, pIDK, pIDS; approx. 5 µg DNA per vial (in buffer solution)

keep at 4°C for short-term storage and in a freezer at -20°C or lower for medium- and long-term storage (take care to avoid repeated freeze-thaw cycles, e.g. by aliquotting DNA prior to freezing)

- **E. coli strains as agar stabs**

a) *E. coli* strain harboring DH10VLP Factory™ bacmid (1 vial)

The DH10VLP Factory™ backbone contains a constitutively expressing mCherry expression cassette which allows for easy monitoring of viral titres via fluorescence without plaque assays.

c) pirHC, pirLC cells<sup>†</sup> (1 vial each)

For propagation and amplification of donor vectors, donor multigene expression constructs or donor-donor fusions

Keep agar stabs at 4°C or at RT; **do not freeze!** We recommend to immediately prepare glycerol stocks from streaked bacterial colonies (see p. 36).

<sup>†</sup> *E. coli* strains expressing the *pir* gene for propagation of donor vectors (any other strain with *pir*<sup>+</sup> background can be used as well). LC: low copy number propagation, HC: high copy number propagation of plasmids with R6Kγ origin.

## Reagents to be supplied by user (see also section D. Protocols)

- Restriction enzymes
- Homing endonucleases PI-SceI and I-CeuI
- Insect cells, e.g. Sf9, Sf21 or High-Five™
- T4 DNA ligase
- Cre recombinase
- Standard *E. coli* strains for cloning (such as TOP10, DH5 $\alpha$ , HB101 etc.)
- Standard laboratory buffers, solutions, media and equipment for bacterial and insect cell culture, transformation etc.
- Commercially available transfection reagents, e.g. FuGENE® (Roche), jetPEI™ (Polyplus transfection), GeneJuice, etc.; alternatively, equipment for electroporation
- Antibiotics, chemicals (e.g IPTG, X-Gal, etc.)

## B. Overview

Virus Like Particles (VLPs) are multiprotein structures that mimic the organization and conformation of authentic native viruses but lack the viral genome, making them non-infectious and safe to handle. VLPs are becoming an increasingly widespread tool for the development of vaccines, and are finding other important applications in biotechnology including serving as a display system for integral membrane proteins for antibody screening, production of immunogens, and for ligand binding assays. The expression of viral structural proteins, such as matrix proteins, can result in the self-assembly of VLPs.

VLP Factory™ is a baculovirus expression system which **constitutively secretes promiscuous influenza virus matrix proteins**, where the system in addition has built in functionality to allow you to introduce either single or multiple additional expression cassettes for production of multiprotein VLPs displaying your surface protein(s), or integral membrane protein(s), of choice. Influenza is an enveloped virus, and VLP Factory™ therefore is a system that constitutively secretes enveloped virus like particles, which will pick up your membrane protein(s) of interest from the insect cell plasma membrane as the particle buds off from the infected cells. Please see **Figure 1** below for a full list of technology applications.

VLP Factory™ transfer vectors are fully synthetic, comprising only functional DNA, and as a result are small and easy to handle, and specifically enable multigene applications (Trowitzsch et al., 2010; Vijayachandran et al, 2011). The VLP Factory™ baculovirus genome (DH10VLP Factory™) is under continuous development to remove harmful/non-essential genes (e.g. v-cath and chiA) and DNA recombination hotspots. This manual also presents a simple and rapid transposition method for integrating your gene(s) of interest into the baculoviral genome and allow for secretion of your protein(s) of interest together with constitutively expressing virus like particles.

Currently, recombinant expression methods require a disproportionate investment in both labor and materials prior to multi-protein expression, and, once expression has been established, provide little or no flexibility for rapidly altering the multiprotein components, which is a prerequisite for revising expression studies. We believe the **VLP Factory™** system will open up new possibilities for all your Virus Like Particle R&D wishes.



**Figure 1.** Applications for VLP produced using VLP Factory™ Technology (colored boxes, top). Electron micrograph of VLP Factory™-produced Tetravalent Flu VLP Containing HA, NA, M1, M2 (bottom left). Electron micrograph of live Flu virus (bottom right).


The baculoviral expression system (BEVS) in its various shapes and forms has become one of the leading eukaryotic expression systems. Baculovirus has become a popular system for a number of applications including the generation of virus-like particles, bioengineering and tissues therapy



research, transduction of mammalian cells, etc. The **VLP Factory™** baculovirus expression system introduced here boasts four major advances that are instrumental in fully exploiting the potential of this heterologous protein production system:


1) The first system that allows you to build your expression needs on top of a constitutive VLP secretion platform. The VLP Factory™ virus constitutively produces a promiscuous influenza VLP which incorporates a wide range of membrane proteins into its envelope. Therefore, if your gene of interest is e.g. a GPCR, expression of your GPCR with the VLP Factory™ will allow you to collect virus like particles from the insect cell culture supernatant displaying your GPCR on the VLP surface.

2) New transfer vectors (pACEBac1, pACEBac2, pIDC, pIDK, pIDS; see Figure 2) that contain a homing endonuclease-based multiplication module. These vectors greatly facilitate modular combination of heterologous genes (in their respective gene expression cassettes) with a minimum requirement for unique restriction sites (BstXI). Baculoviral promoters (currently p10 and polh very late promoters) can be exchanged in our vectors for other promoter sequences (early and late viral, mammalian) if desired. Likewise, terminator sequences (currently SV40, HSVtk) can be substituted as required.

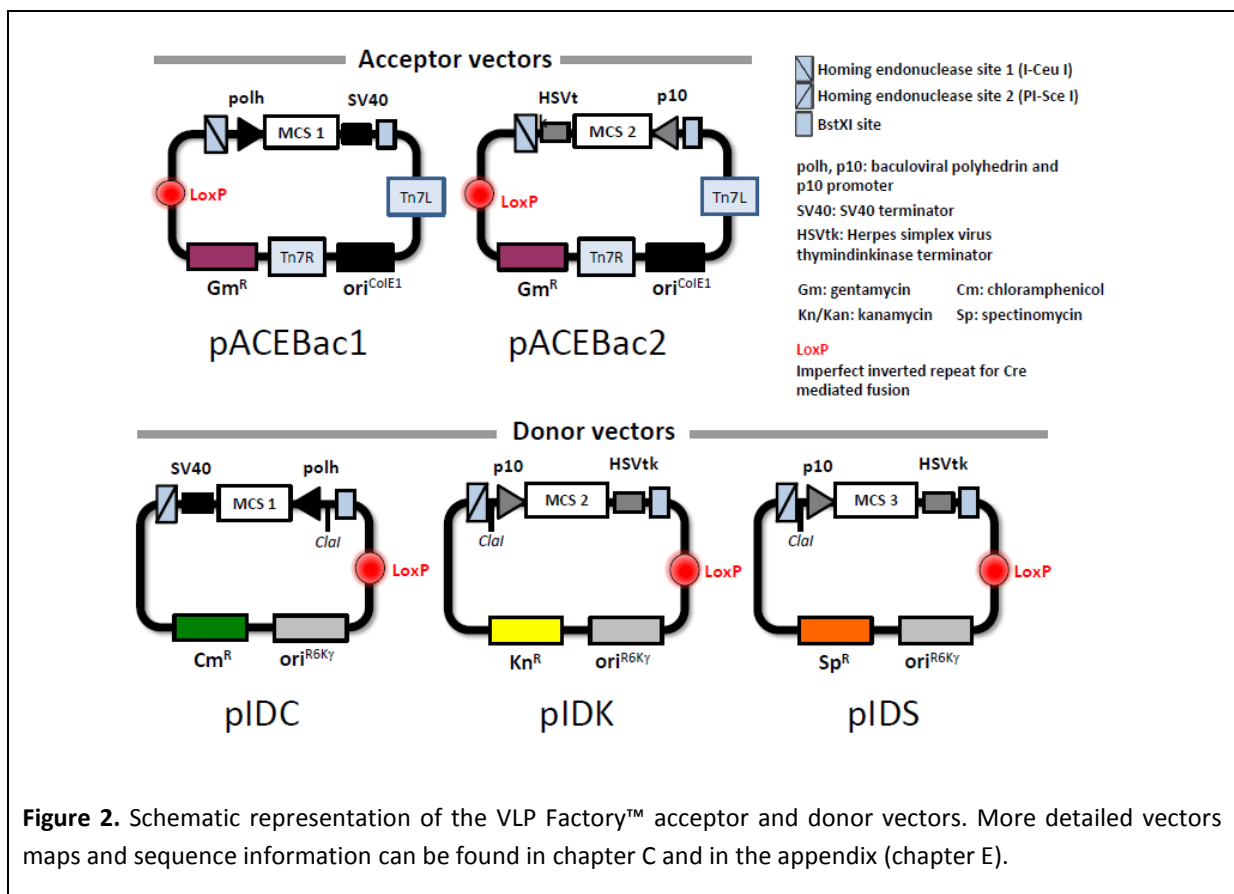


**3)** The baculovirus genome has been engineered to improve its protein production properties. Disruption of two baculoviral genes improves the integrity of cellular compartments during infection and protein production (see below). The quality of proteins produced by this system is significantly improved by curtailing virus-dependent proteolytic activity and reducing cell lysis.

**4)** New protocol for rapid generation of multigene expression constructs via Cre-LoxP recombineering. The resulting multigene fusion is then incorporated into baculovirus DNA by accessing the viral genome via a specific site (see Figure 3). This protocol can be used to integrate multigene cassettes with coding sequences for multiprotein complex subunits into VLP Factory™, but also to integrate specific enzymes (kinases, acetylases etc.) for modifying the proteins under investigation.

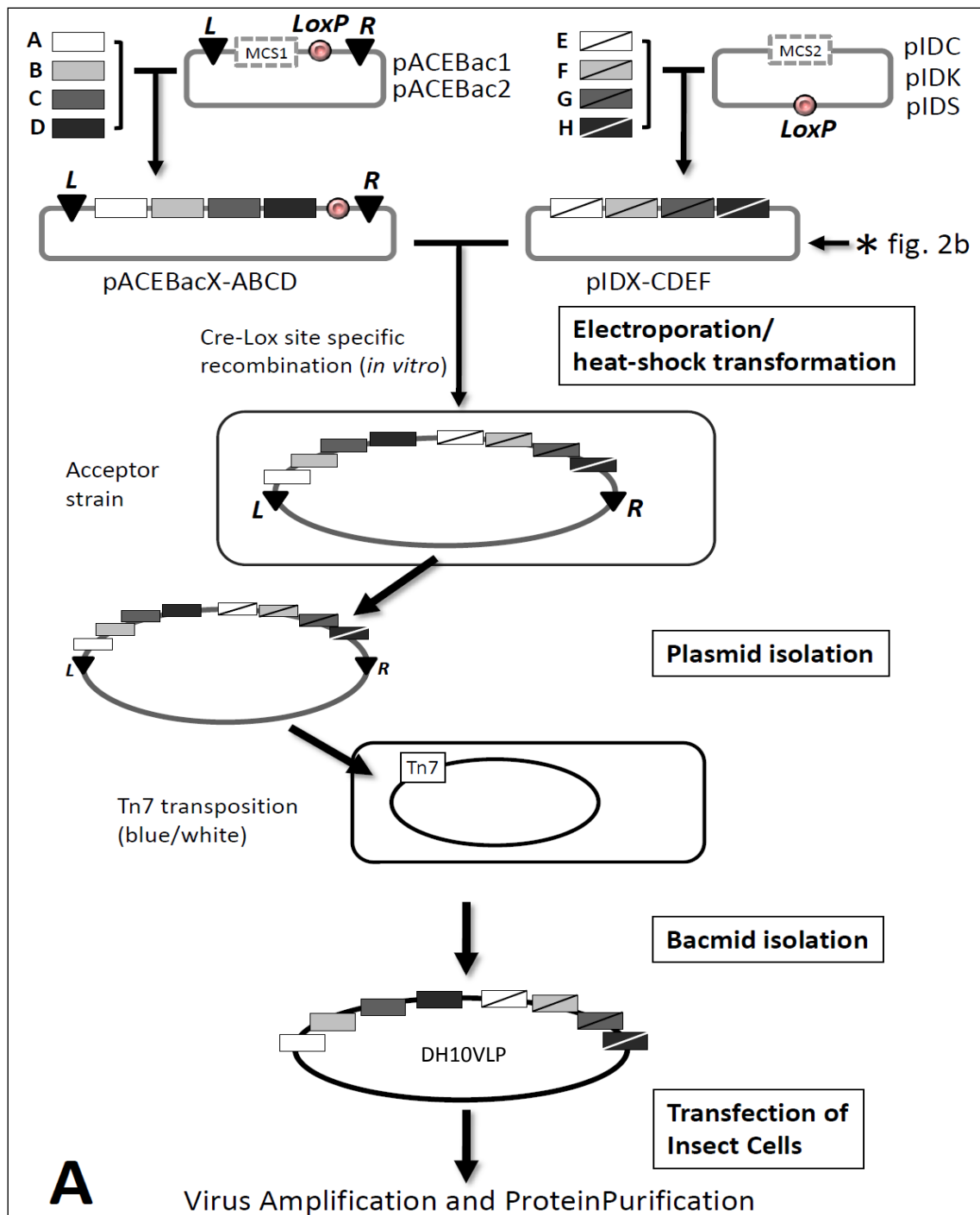

**Figure 3. Schematic overview of the VLP Factory™ system and its application.** Genes of interest are assembled into multigene expression cassettes using the multiplication module present on the donor (pIDC, pIDK, pIDS) and acceptor vectors (pACEBac1, pACEBac2). Acceptor-donor fusions can then be generated by Cre-LoxP recombination. These multigene fusions contain one Acceptor and one to several Donor vectors, each with one or several genes of interest (here A-H). Desired Acceptor-Donor combinations are identified by transforming in E.coli and subsequent antibiotic selection. Acceptor vectors contain the DNA elements required for integration into the baculovirus via Tn7 transposition. DH10VLP Factory™ E. coli cells contain the recipient baculovirus and also the transposase required for Tn7 transposition for integrating acceptor and acceptor-

donor derivatives that are transformed into these cells. Colonies containing bacmid carrying integrated multigenecassettes are identified by blue/white screening (Tn7 transposition disrupts expression of the lacZ $\alpha$  peptide) plus gentamycin resistance, and optionally through selective resistance markers introduced via Cre-catalyzed integration and hosted on the acceptor, donor or acceptor-donor derivatives). Bacmid DNA is prepared from selected clones and used to transfect insect cells for VLP protein production. LoxP sites in the acceptor-donor fusions have been omitted for reasons of clarity.

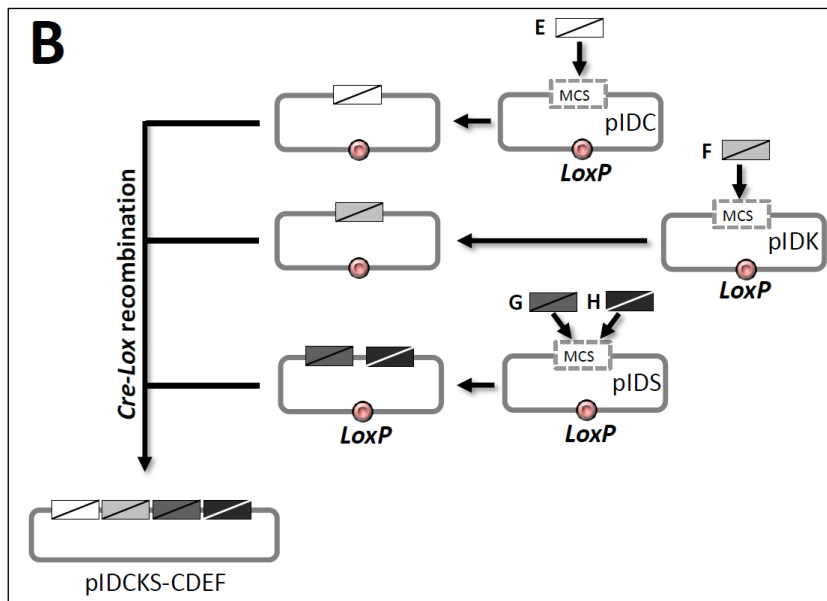


**Figure 4. Generation of multigenedonor constructs through Cre-Lox fusion.** As indicated in Figure 3, donor multigeneexpression cassette constructs can also be generated by Cre-Lox recombination. Individual or multiple gene cassettes are cloned into the multiple cloning site via standard restriction-ligation cloning or, when introducing multiple gene cassettes, homing endonuclease /BstXI cloning. The gene cassettes harbored on different donor vectors are then merged into a single vector construct via Cre-Lox recombination. This construct will differ from the multigeneconstructs in Figure 2a with respect to selective markers. While the multigeneconstruct in Figure 3 carries only one antibiotic resistance marker, the construct here in Figure 4 will carry three, one from each donor vector. This will allow selection of multigeneconstructs with higher stringency by subjecting the constructs to a multi-antibiotic selection regimen (refer to protocol 2). LoxP sites in the donor fusion have been omitted for reasons of clarity.

## C. New Baculovirus Tools for for Virus Like Particle (VLP) Applications

### C.1. Transfer vectors: the Acceptor-Donor recombineering system.

The **Acceptor vectors** pACEBac1 and pACEBac 2 contain multiple cloning sites (see Appendix) flanked by either a polh or p10 promoters and SV40 or HSVtk polyA signal sequences, respectively. A multiplication module M – defined by the homing endonuclease site I-CeuI and a corresponding BsXI site (see Figure 5) – allows integration of multiple gene cassettes (ORFs and associated regulatory regions). The sequences used for Tn7 transposition (Tn7L and Tn7R) encompass the expression cassettes and a gentamycin resistance marker.

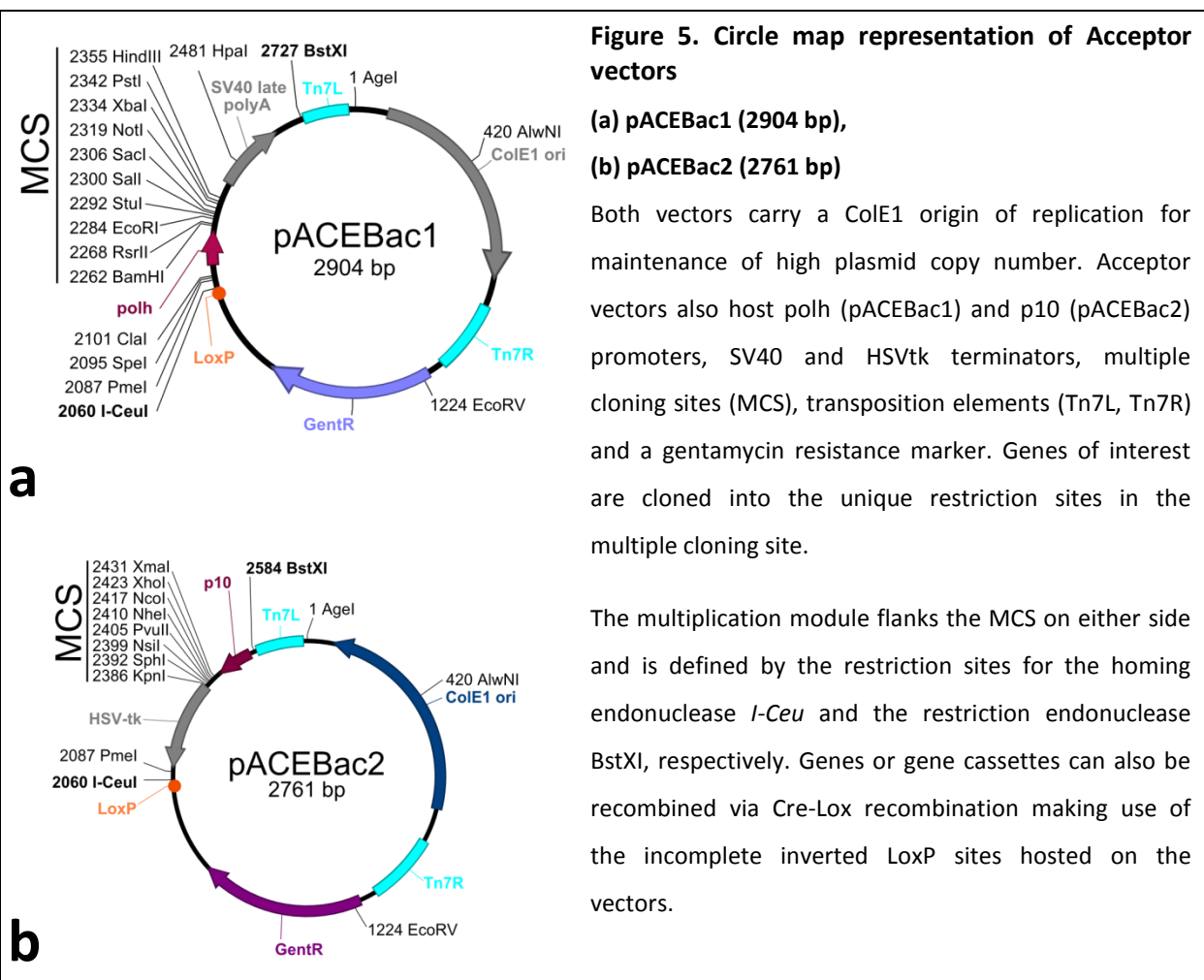


The **Donor vectors** pIDC, pIDS, pIDK are similar to the acceptor vectors with respect to their over-all design. The multiple cloning site is bracketed by a multiplication element (in this case, PI-SceI /BstXI) to enable concatenation of inserts between the different donor vectors. Vectors also contain a LoxP incomplete inverted repeat to create acceptor-donor or donor-donor fusions. The vectors contain “tell-tale” resistance markers (pIDC: chloramphenicol, pIDK: kanamycin, pIDS: spectinomycin) and, importantly, a conditional R6Ky origin of replication which makes its propagation dependent on the expression of the *pir* gene in the prokaryotic host (such as the pirLC and pirHC cells contained in the kit).

**Figure 6. Circle map representation of Donor vectors**

**a) pIDC, b) pIDK, c) pIDS.**

Circle maps show promoters (polh, p10), terminators (SV40, HSVtk), multiple cloning sites (MCS), the incomplete inverted repeat for *cre-lox* site-specific recombination (LoxP) and a resistance marker (chloramphenicol, kanamycin, and spectinomycin, respectively). Genes of interest are cloned into the MCS using unique restriction sites. The multiplication module consists of the homing endonuclease site PI-SceI and the restriction endonuclease site BstXI.

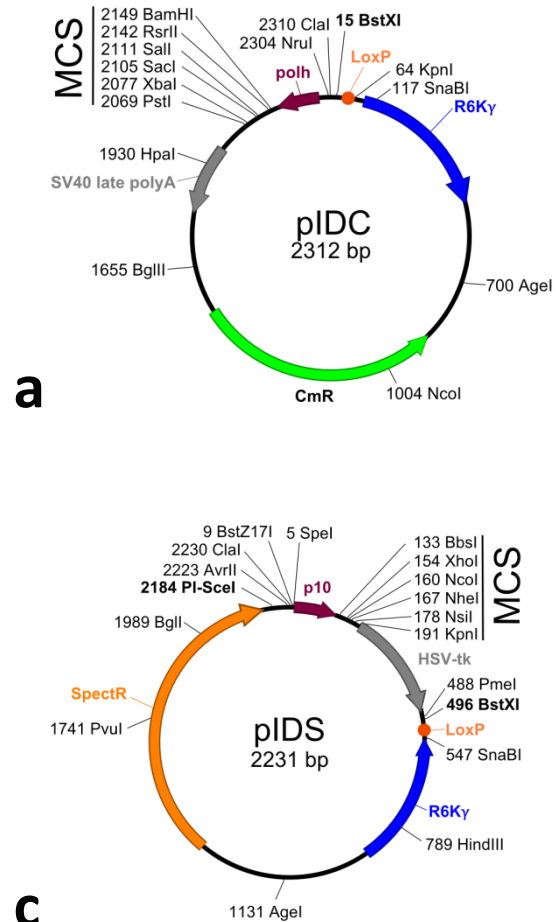

Currently, the VLP Factory™ vectors do not contain DNA sequences coding for affinity tags that will facilitate purification or solubilization of the protein(s) of interest. Tags that are typically used are C- or N-terminal oligohistidine tags, with or without protease sites for tag removal. They can be introduced by designing the respective PCR primers used for amplification of the genes of interest. We recommend outfitting Donors or Acceptors of choice with any custom tag that is favored in individual

user laboratories prior to inserting recombinant genes of interest. This is best done by using a design that will, after tag insertion, still be compatible with the recombination-based principles of ACEMBL system usage.

## C.2. Generating multigene expression cassettes

### C.2.1. Multigene construction via homing endonuclease/*Bst*XI multiplication


The acceptor and donor vectors are suited for generating multigene expression cassettes from individual gene expression cassettes (complete with regulatory regions such as promoter and terminator) via a multiplication module bracketing the multiple cloning site (MCS). All VLP Factory™ vectors contain a homing endonuclease (HE) site and a matching designed *Bst*XI site that envelop the MCS. Homing endonucleases have long recognition sites (20-30 base pairs or more). Although not all equally stringent, homing endonuclease sites are most probably unique in the context of even large plasmids, or, in fact, entire genomes.

The logic of multiplication is illustrated below. The homing endonuclease site can be used to insert entire expression cassettes into a vector already containing one gene or several genes of interest as separate expression cassettes. The only prerequisite for assembling multigene expression cassettes is that the homing endonucleases and restriction enzymes used for multiplication (*I-CeuI*/*PI-SceI* and *Bst*XI) are unique, which can be easily accomplished for instance by site directed mutagenesis prior to multigene cassette assembly. First, individual genes are cloned into the multiple cloning sites of the acceptor and donor vectors. The entire expression cassette, including promoter and terminator, is then excised by *I-CeuI* and *Bst*XI (acceptors) or *PI-SceI* and *Bst*XI (donors) digestion. The resulting fragment is placed into the multiplication module of another acceptor or donor vector containing single or multiple gene cassettes. The restriction sites involved are eliminated in the process and multiplication can be repeated iteratively using the module present in the inserted cassette. Moreover, promoter and terminator sequences can be easily modified if desired using appropriate restriction sites in our vectors.




**Please note that multiplication cannot be accomplished from donors to acceptors and vice versa since the overhangs generated by endonuclease digestion are incompatible.**



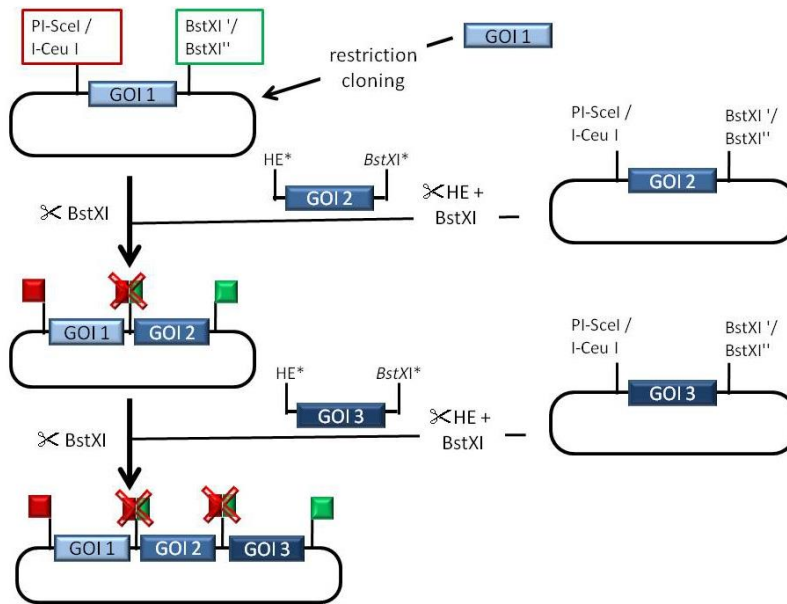

**Figure 7. Assembling individual gene cassettes into multigene expression cassettes.** The logic of multiplication is shown schematically. The expression cassette containing the gene of choice (denoted as GOI2 in this case) is excised by digestion with the homing endonuclease (red box) and BstXI (green box). For acceptors vectors, I-CeuI is the homing endonuclease of choice, and for donor vectors PI-SceI. The plasmid vector harboring the GOI1-cassette only needs to be linearized with BstXI. The homing endonucleases produce cohesive ends that are compatible with the ends generated by the BstXI digest. Upon insertion of GOI2 into the target vector, a homing endonuclease/BstXI hybrid restriction site is created that can then cannot be cut by either enzyme (crossed-out red/green box) while the 3'-BstXI site is regenerated. The same procedure can be repeated over and over as exemplified by the integration of GOI3. This cycling logic can be used to generate multigeneassemblies. Note that the promoters and terminators are not explicitly shown for reasons of clarity.

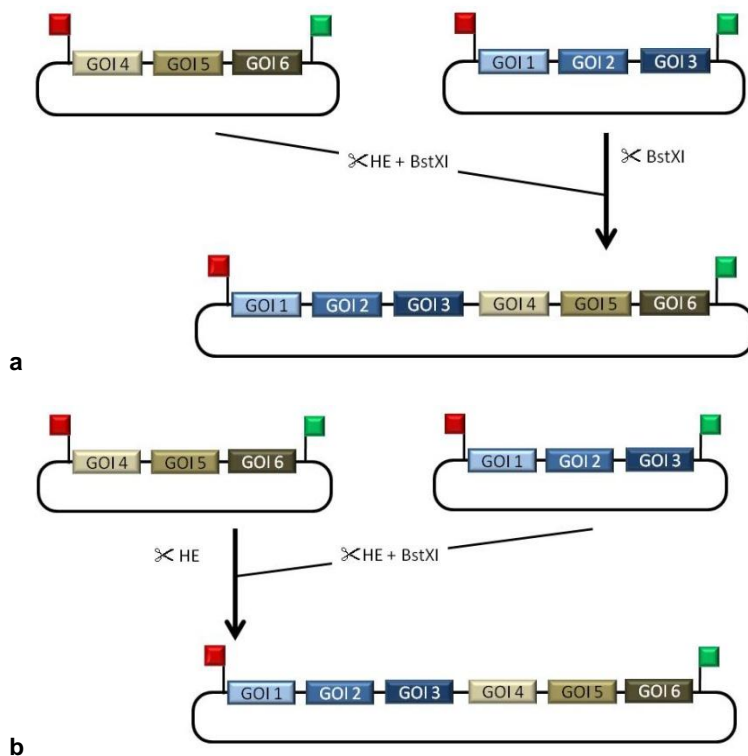

**Figure 8. Combining multigene expression cassettes.** Different multigeneexpression cassettes can be combined into one expression construct following the same logic that applies to the generation of multigeneexpression cassettes from individual gene cassettes. The 5' homing endonuclease recognition site (filled red box) will be preserved if GOI1 has been introduced by conventional restriction cloning into the MCS. Promoters and terminators are not explicitly shown for reasons of clarity but flank the GOIs in every individual gene expression cassette.

### C.2.2. Multigene construction using Cre-Lox recombination

Cre recombinase is a member of the integrase family (Type I topoisomerase from bacteriophage P1). It recombines a 34 bp loxP site in the absence of accessory protein or auxiliary DNA sequence. The loxP site is comprised of two 13 bp recombinase-binding elements arranged as inverted repeats which flank an 8 bp central region where cleavage and the ligation reaction occur.


The site-specific recombination mediated by Cre recombinase involves the formation of a Holliday junction (HJ). The recombination events catalyzed by Cre recombinase depend on the location and relative orientation of the loxP sites. Two DNA molecules, for example an acceptor and a donor plasmid, containing single loxP sites will be fused. Furthermore, the Cre recombination is an equilibrium reaction with 20-30% efficiency in recombination. This provides useful options for multigenecombinations for multi-protein complex expression.

13bp	8bp	13bp
5'...ATAACTTCGTATA	GCATACAT	TATACGAAGTTAT...3'
3'...TATTGAAGCATAT	CGTATGTA	ATATGCTTCAATA...5'
inverted repeat	Spacer	inverted repeat

**Figure 9.: LoxP imperfect inverted repeat**

In a reaction where several DNA molecules such as donors and acceptors are incubated with Cre recombinase, the fusion/excision activity of the enzyme will result in an equilibrium state where single vectors (educt vectors) and all possible fusions coexist. Donor vectors can be used with acceptors and/or donors, and vice versa. Higher order fusions are also generated where more than two vectors are fused. This is shown schematically in Figure 10 below.

The fact that Donors contain a conditional origin of replication that depends on a *pir*<sup>+</sup> (*pir* positive) background now allows for selecting out from this reaction mix all desired Acceptor-Donor(s) combinations. For this, the reaction mix is used to transform *pir* negative strains (TOP10, DH5 $\alpha$ , HB101 or other common laboratory cloning strains). Then, Donor vectors will act as suicide vectors when plated out on agar containing the antibiotic corresponding to the Donor encoded resistance marker, unless fused with an Acceptor. By using agar with the appropriate combinations of antibiotics, all desired Acceptor-Donor fusions can be selected for.

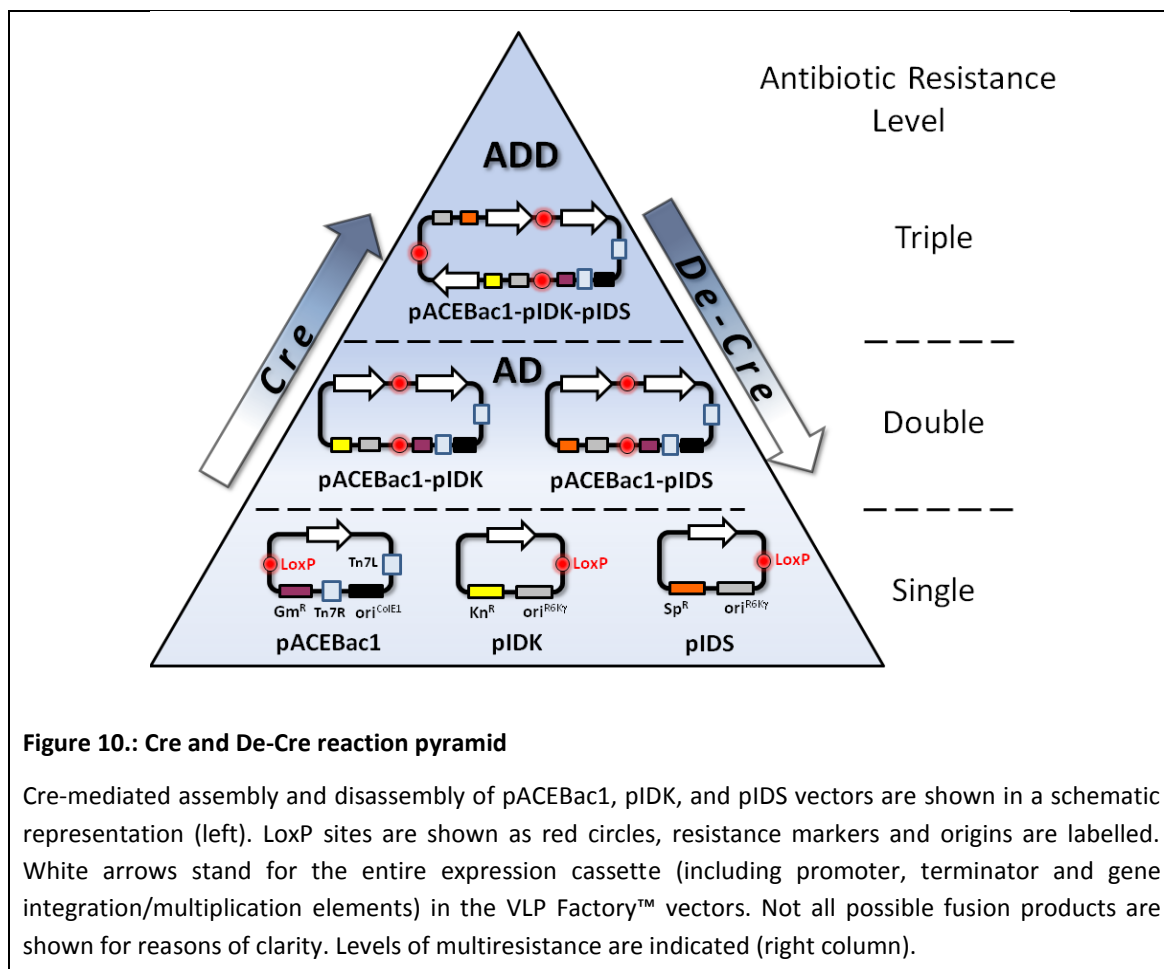


### C.2.3. Combining HE/BstXI cycling and Cre-Lox recombination

Of course, both methods can also be combined to generate multiple gene-expression cassette constructs. To this end, you can introduce multiple gene cassettes with the homing endonuclease/BstXI protocol into different Acceptor/Donor vectors and then fuse these using the Cre-Lox modules.

### C.3. Baculovirus engineered for improved protein production.

During heterologous protein production using other commercially available baculovirus expression system, viral-dependent proteolytic breakdown consistent with the action of a cysteine protease can be observed. The VLP Factory™ baculovirus genome (schematically shown in Figure 11) was modified to yield improved protein production properties. Two baculoviral genes, *v-cath* and *chiA*, have been disrupted which leads to improved maintenance of cellular compartments during infection and protein production. The *v-cath* gene (Slack et al., 1995) encodes for a viral cathepsin-type cysteine protease, V-CATH, which is activated upon cell death by a process that depends on a juxtaposed gene on the viral DNA, *chiA*, which encodes for a chitinase (Hawtin et al., 1995; Hom and Volkman, 2000). Both are involved in the liquefaction of the host insect cells (Slack et al., 1995; Hawtin et al., 1997).

## VLP Factory™ Baculoviral D

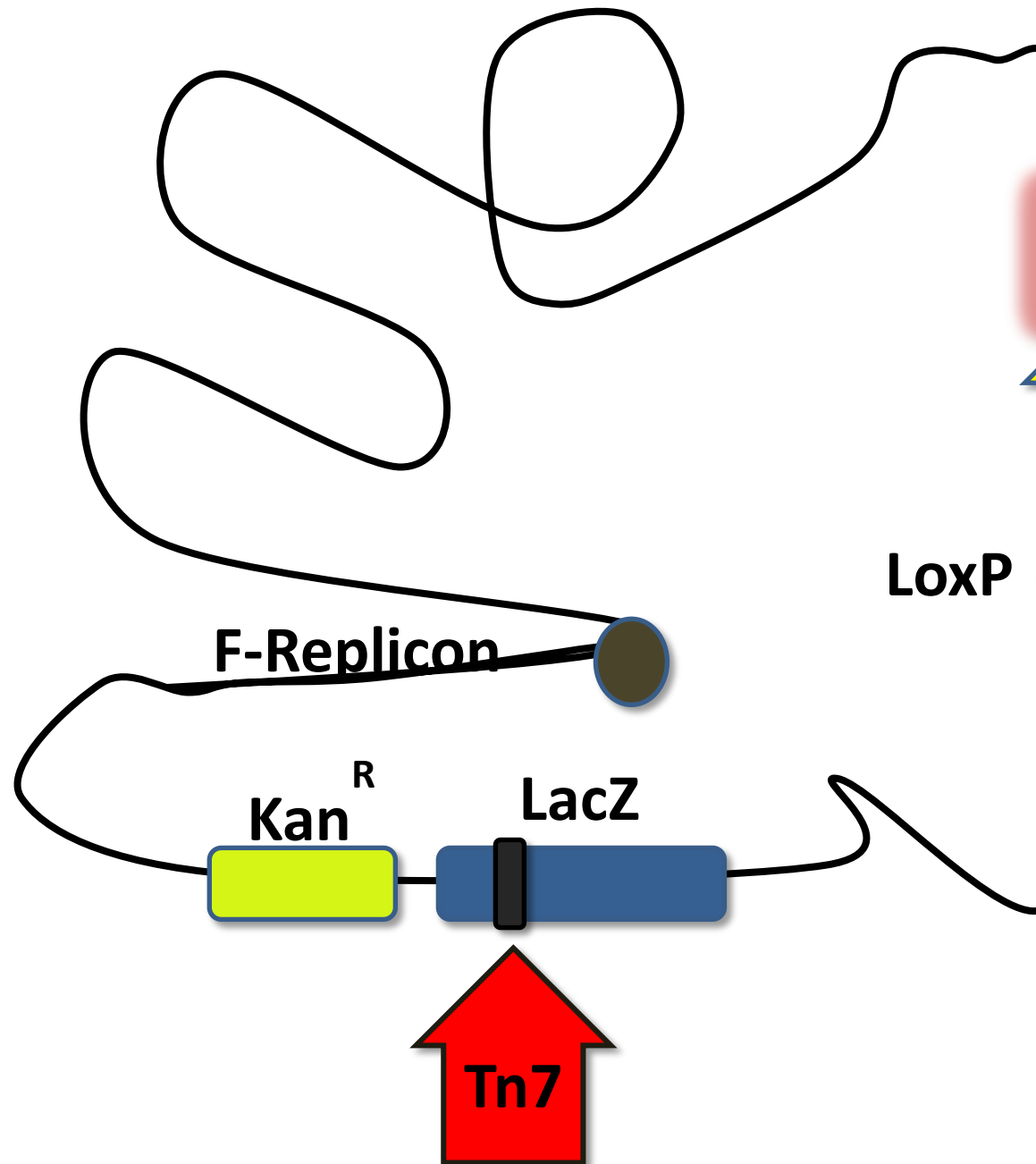

- ✓ Co-expression of Flu M
- ✓ Specific Viral Gene Del
- ✓ Two Separate Viral Ent
- ✓ Virally Encoded Fluore

**Figure 11.: DH10VLP Factory™ baculoviral DNA.** The modified viral genome is shown in a schematic representation. The Tn7 attachment site is located within a *LacZα* gene; insertion of Tn7 elements from pACEBac derivatives therefore produces a white phenotype when plated on agar containing BluoGal and IPTG. The viral genes *v-cath* and *chiA* are disrupted by replacement with an antibiotic marker and a non-functional LoxP sequence (both not shown for reasons of clarity).

Disruption of both genes served to a) eliminate V-CATH activity and b) to enable chitin-affinity chromatography for purification without interference from the *chiA* gene product. The quality of proteins produced by the VLP Factory™ baculovirus is significantly improved through a reduction of virus-dependent proteolytic activity and reduced cell lysis. The disrupted viral DNA sequence was replaced with a LoxP sequence for *cre-lox* site-specific recombination. Note that this LoxP site is not used for introduction of the target constructs into VLP Factory™. Instead, the genes of interest are transferred to the bacmid via transposition into the mini Tn7 attachment site. Successful integration results in disruption of the *lacZα* subunit-coding sequence. As a consequence, clones carrying inserted DNA will appear white.

## C.4. Introducing additional control elements

Protein expression requires certain genetic elements at the DNA and RNA level for it to work efficiently, as exemplified by the Shine-Dalgarno or Kozak sequence in bacterial and vertebrate mRNAs, respectively. Without this ribosome-binding motif, efficient translation of the corresponding mRNAs is compromised. Yet, many genes/mRNAs lack such a consensus sequence motif which is not an absolute must for successful translation.

While the polyhedrin promoter already is one of the strongest promoters known, protein production may in some cases be enhanced by substituting the intrinsic leader with alternative leader sequences in the gene expression cassettes. While there is no guarantee that this will work equally well for all proteins under investigation, several reports suggest the inclusion of such motifs into gene design, e.g. the L21 leader from a lobster tropomyosin cDNA (Sano et al., 2002).

In most cases, though, protein production should work just fine with the promoters provided in the transfer vectors.

## D. Protocols

### D.0 Introductory remarks

Please note that the bacteria in the agar stabs have not been made competent for transformation. If you wish to use them to transform your constructs, you will have to prepare competent cells. This applies specifically to the pirHC and pirLC strains used to maintain and amplify donor constructs. You may follow your preferred protocol for preparing chemically or electrocompetent cells, e.g. Inoue et al. (1990) or variations of this protocol, or standard protocols as described in Current Protocols in Molecular Biology or Sambrook and Russell: Molecular Cloning (3<sup>rd</sup> edition, 2001, or older versions).

### D.1 Cloning into pACEBac or pIDx transfer vectors

Reagents:

Restriction endonucleases

DNA ligase

E. coli competent cells

Antibiotics: Chloramphenicol, Gentamycin, Kanamycin, Spectinomycin

The genes of choice are cloned using standard cloning procedures into the multiple cloning sites MCS (see *Supplementary Information*) of pACEBac1/2 and pIDC, pIDK, pIDS. Ligation reactions for pACEBac derivatives are transformed into standard *E. coli* cells for cloning (such as TOP10, DH5 $\alpha$ , HB101) and plated on agar containing gentamycin (7  $\mu$ g/ml). Ligation reactions for pIDx derivatives are transformed into *E. coli* cells expressing the *pir* gene (pirHC and pirLC from this kit – in this case you will need to make the cells electro- or chemically competent first; other strains, e.g. BW23473, BW23474) and plated on agar containing chloramphenicol (pIDC; 25  $\mu$ g/ml), kanamycin (pIDK; 50  $\mu$ g/ml) or spectinomycin (pIDS; 50  $\mu$ g/ml). Correct clones are selected based on specific restriction digestion patterns and DNA sequencing of the inserts.

### D.2 Multiplication by using the HE and BstXI sites

VLP Factory™ donor vectors contain a recognition site for the homing endonuclease PI-SceI (Figures 5-6). Upon cleavage, this HE site yields a 3' overhang with the sequence -GTGC. Acceptor vectors contain



the homing endonuclease site I-CeuI (see Figure 5-6), which upon cleavage will result in a 3' overhang of -CTAA. On acceptors and donors, the respective HE site precedes the MCS (see Figures 7-8). The 3' end of the MIE contains a specifically designed BstXI site, which upon cleavage will generate a matching overhang. The basis of this is the specificity of cleavage by BstXI. The recognition sequence of BstXI is defined as CCANNNNN'NTGG (the apostrophe marks the position of the phosphodiester link cleavage). The residues denoted as N can be chosen freely. Donor vectors thus contain a BstXI recognition site with the sequence CCATGTGC'CTGG, and Acceptor vectors contain CCATCTAA'TTGG. The overhangs generated by BstXI cleavage in each case will match the overhangs generated by HE cleavage. Note that Acceptors and Donors have different HE sites.

The recognition sites are not symmetric. Therefore, ligation of a HE/BstXI digested fragment into a HE site of an VLP Factory™ vector will be (1) directional and (2) result in a hybrid DNA sequence where a HE halfsite is combined with a BstXI half site (see Figure 5). This site will be cut by neither the HE nor BstXI. Therefore, in a construct that has been digested with a HE, insertion by ligation of HE/BstXI digested DNA fragment containing an expression cassette with one or several genes will result in a construct which contains all heterologous genes of interest, enveloped by an intact HE site in front, and a BstXI site at the end. Therefore, the process of integrating entire expression cassettes by means of HE/BstXI digestion and ligation into a HE site can be repeated iteratively.

## **D2.1 Protocol 1. Multiplication by using homing endonuclease/BstXI.**

### Reagents required:

Homing endonucleases PI-SceI, I-CeuI  
10x Buffers for homing endonucleases  
Restriction enzyme BstXI (and 10x Buffer)  
T4 DNA ligase (and 10x Buffer)  
*E. coli* competent cells  
Antibiotics

### **Step 1: Insert preparation**

Restriction reactions are carried out in 40 µl reaction volumes, using homing endonucleases PI-SceI (Donors) or I-CeuI (Acceptors) as recommended by the supplier.

Acceptor or donor plasmid ( $\geq 0.5 \mu\text{g}$ ) in ddH <sub>2</sub> O	32 $\mu\text{l}$	
10x restriction enzyme buffer	4 $\mu\text{l}$	
10 mM BSA	2 $\mu\text{l}$	
PI-SceI (Donors) or I-CeuI (acceptors)	2 $\mu\text{l}$	ca. 10 U

Reactions are then purified using a PCR extraction kit or by acidic ethanol precipitation, and subsequently digested with BstXI according to the supplier's recommendations.

HE digested DNA in ddH <sub>2</sub> O	32 $\mu\text{l}$	
10x restriction enzyme buffer	4 $\mu\text{l}$	
10 mM BSA	2 $\mu\text{l}$	
BstXI	2 $\mu\text{l}$	ca. 20 U

Gel extraction of insert(s):

Processed insert is then purified by agarose gel extraction using commercial kits (Qiagen, MachereyNagel etc). Elution of the extracted DNA in the minimal volume defined by the manufacturer is recommended.

## Step 2: Vector preparation

Restriction reactions are carried out in 40  $\mu\text{l}$  reaction volumes, using homing endonucleases PI-SceI (Donors) or I-CeuI (Acceptors) as recommended by the supplier.

Acceptor or donor plasmid ( $\geq 0.5 \mu\text{g}$ ) in ddH <sub>2</sub> O	33 $\mu\text{l}$	
10x Restriction enzyme buffer	4 $\mu\text{l}$	
10 mM BSA	2 $\mu\text{l}$	
BstXI	2 $\mu\text{l}$	ca. 20 U

Reactions are then purified by PCR extraction kit or acidic ethanol precipitation, and next treated with intestinal alkaline phosphatase according to the supplier's recommendations. Dephosphorylation is performed to minimize vector re-annealing and to increase integration of the insert.

HE digested DNA in ddH <sub>2</sub> O	17 $\mu\text{l}$	
10x Alkaline phosphatase buffer	2 $\mu\text{l}$	
Alkaline phosphatase	1 $\mu\text{l}$	ca. 10 U

Gel extraction of vector:

Processed vector is then purified by agarose gel extraction using commercial kits (Qiagen, MachereyNagel etc). Elution of the extracted DNA in the minimal volume defined by the manufacturer is recommended.

### Step 3: Ligation

Ligation reactions are carried out in 20 µl reaction volumes:

HE/Phosphatase treated vector (gel extracted)	4 µl	
HE/BstXI treated insert (gel extracted)	14 µl	
10x T4 DNA Ligase buffer	2 µl	
T4 DNA Ligase	0.5 µl	150 - 200 c.e. U or 2-4 Weiss U

Ligation reactions are performed at 25°C for 1h or at 16°C overnight.

### Step 4: Transformation

Mixtures are next transformed into competent cells following standard transformation procedures.

Ligation reactions for pACEBac1 and pACEBac2 derivatives are transformed into standard *E. coli* cells for cloning (such as TOP10, DH5α, HB101) and, after recovery, are plated on agar containing gentamycin (7 µg/ml).

Reactions for Donor derivatives are transformed into *E. coli* cells expressing the *pir* gene (such as BW23473, BW23474, or PIR1 and PIR2 from Invitrogen and, of course, pirLC and pirHC in this kit) and plated on agar containing chloramphenicol (25 µg/ml, pIDC), kanamycin (50 µg/ml, pIDK), or spectinomycin (50 µg/ml, pIDS).

### Step 5: Plasmid analysis

Plasmids are cultured and correct clones selected based on specific restriction digestion and DNA sequencing of the inserts.

## D.3 Cre-LoxP reaction of Acceptors and Donors

### D 3.1. Protocol 2: Cre-LoxP fusion of Acceptors and Donors

This protocol is designed for generating multigene fusions from Donors and Acceptors by Cre-LoxP reaction.

Reagents:

Cre recombinase (from NEB or self made)

Standard *E. coli* competent cells (*pir*<sup>-</sup> strain)

Antibiotics

96well microtiter plates

12 well tissue-culture plates (or petri dishes) w. agar/antibiotics

LB medium

1. For a 20µl Cre reaction, mix 1-2 µg of each educt in approximately equal amounts. Add ddH<sub>2</sub>O to adjust the total volume to 16-17 µl, then add 2 µl 10x Cre buffer and 1-2µl Cre recombinase (1-2 U).
2. Incubate Cre reaction at 37°C (or 30°C) for 1 hour.
3. Optional: load 2-5 µl of Cre reaction on an analytical agarose gel for examination.

*Heat inactivation at 70°C for 10 minutes before gel loading is strongly recommended.*

4. For chemical transformation, mix 10-15µl Cre reaction with 200 µl chemically competent cells. Incubate the mixture on ice for 15-30 minutes. Then perform heat shock at 42°C for 45-60 s.

*Up to 20 µl Cre reaction (0.1 volumes of the chemically competent cell suspension) can be directly transformed into 200 µl chemical competent cells.*

For electrotransformation, up to 2 µl Cre reaction can be directly mixed with 100 µl electrocompetent cells, and transformed by using an electroporator (e.g. BIORAD *E. coli* Pulser) at 1.8-2.0 kV.

*Larger volumes of Cre reaction must be desalted by ethanol precipitation or via PCR purification columns before electrotransformation. The desalted Cre reaction mix should not exceed 0.1 volumes of the electrocompetent cell suspension.*

*The cell/DNA mixture can be immediately used for electrotransformation without prolonged incubation on ice.*

5. Add up to 400 µl of LB (or SOC) medium per 100 µl of cell/DNA suspension immediately after the transformation (heat shock or electroporation).

6. Incubate the suspension in a 37°C shaking incubator overnight or for at least 4 hours (recovery period).

*To recover multifusion plasmid containing more than 2 resistance markers, it is strongly recommended to incubate the suspension at 37°C overnight.*

7. Plate out the recovered cell suspension on agar containing the desired combination of antibiotics. Incubate at 37°C overnight.

8. Clones from colonies present after overnight incubation can be verified by restriction digestion at this stage (refer to steps 12-16).

*This quality control step should be carried out especially in the case that only one specific multifusion plasmid is desired.*

For further selection by single antibiotic challenges on a 96 well microtiter plate, continue to step 9.

*Several to many different multifusion plasmid combinations can be processed and selected in parallel on one 96 well microtiter plate.*

9. For 96 well antibiotic tests, inoculate four colonies from each agar plate with different antibiotic combinations into approx. 500 µl LB medium without antibiotics. Incubate the cell cultures in a 37°C shaking incubator for 1-2 hours.


10. While incubating the colonies, fill a 96 well microtiter plate with 150 µl antibiotic-containing LB medium. It is recommended to add coloured dye (positional marker) in the wells indicated.

*A typical arrangement of the solutions, which is used for parallel selections of multifusion plasmids, is shown in Figure 12. The concept behind the 96 well plate experiment is that every cell suspension from single colonies needs to be challenged by all four single antibiotics for unambiguous interpretation.*

11. Add 1 µl aliquots of pre-incubated cell culture (Step 9) to the corresponding wells. Then incubate the inoculated 96 well microtiter plate in a 37°C shaking incubator overnight at 180-200 rpm.

*Recommended: use parafilm to wrap the plate to avoid drying out.*

*The remainder of the pre-incubated cell cultures can be kept at 4°C for further inoculations if necessary.*

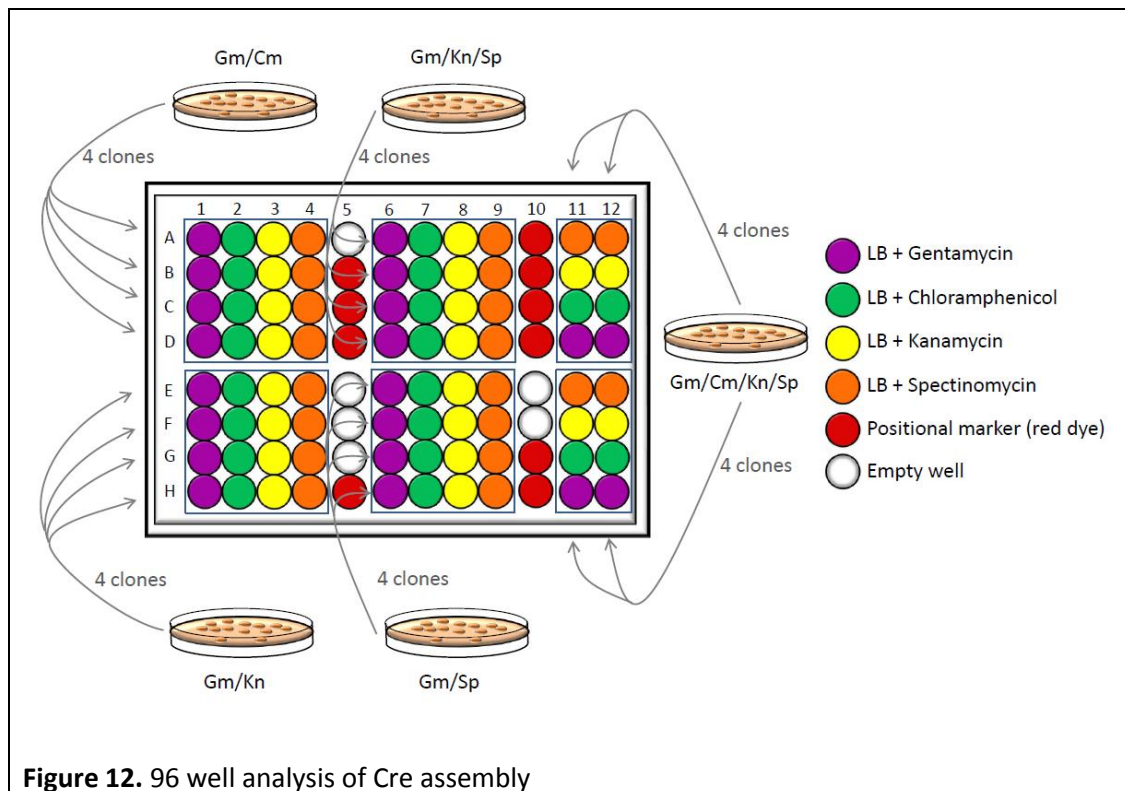


**Figure 12.** 96 well analysis of Cre assembly

12. Select transformants containing desired multifusion plasmids based on antibiotic resistance, according to the combination of dense (positive) and clear (no growth) cell microcultures from each colony. Inoculate 10-20  $\mu$ l cell culture into 10 ml LB media with corresponding antibiotics. Incubate in a 37°C shaking incubator overnight.
13. Centrifuge the overnight cell cultures at 4000g for 5-10 minutes. Purify plasmid from the resulting cell pellets with common plasmid miniprep kits, according to manufacturers' recommendation.
14. Determine the concentrations of purified plasmid solutions by using UV absorption spectroscopy (e.g. by using a NanoDrop™ 1000 machine).
15. Digest 0.5-1  $\mu$ g of the purified plasmid solution in a 20  $\mu$ l restriction digestion with appropriate endonuclease(s). Incubate under recommended reaction condition for approx. 2 hours.
16. Use 5-10  $\mu$ l of the digestion for analytical agarose (0.8-1.2%) gel electrophoresis. Verify plasmid integrity by comparing the experimental restriction pattern to a restriction pattern predicted *in silico* (e.g. by using program VectorNTI from Invitrogen or similar programs).

### D 3.2. Protocol 3. Deconstruction of fusion vectors by Cre

The following protocol is suitable for releasing any single educt from multifusion constructs (deconstruction). This is achieved by Cre-LoxP reaction, transformation and plating on agar with appropriately reduced antibiotic resistance level (cf. Figure 10). In the liberated educt entity, encoding genes can be modified and diversified. Then, the diversified construct is resupplied by Cre-LoxP reaction (C3.1).

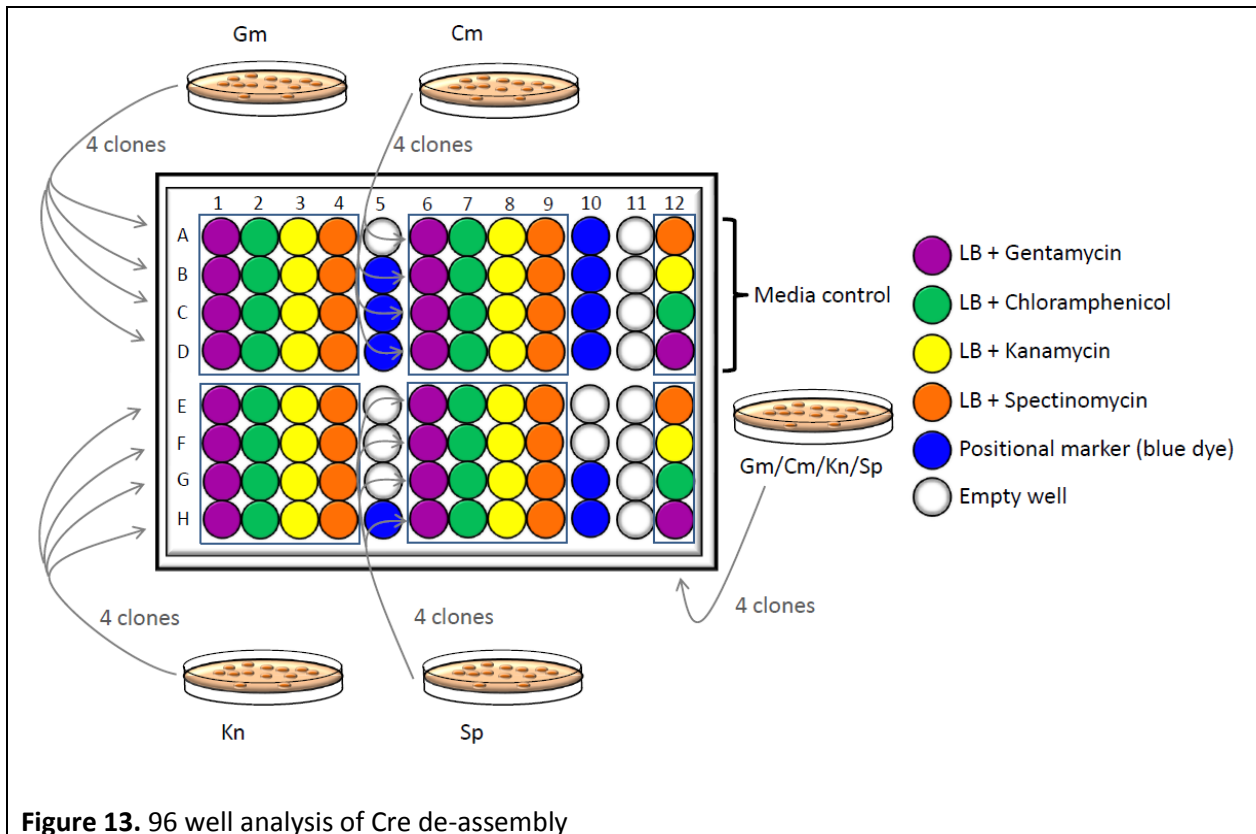

**Figure 13.** 96 well analysis of Cre de-assembly

#### Reagents:

Cre recombinase (and 10x Buffer)

*E. coli* competent cells

(*pir*<sup>+</sup> strains, *pir*<sup>-</sup> strains can be used only when partially deconstructed Acceptor-Donor fusions are desired).

Antibiotics

1. Incubate approx. 1 µg multifusion plasmid with 2 µl 10x Cre buffer, 1-2 µl Cre recombinase, add ddH<sub>2</sub>O to adjust the total reaction volume to 20 µl.
2. Incubate this Cre deconstruction reaction mixture at 30°C for 1 to 4 hour(s).

3. Optional: load 2-5 µl of the reaction on an analytical agarose gel for examination.

*Heat inactivation at 70°C for 10 minutes before gel loading is strongly recommended.*

4. For chemical transformation, mix 10-15µl De-Cre reaction with 200 µl chemically competent cells. Incubate the mixture on ice for 15-30 minutes. Then perform heat shock at 42°C for 45-60 s.

*Up to 20 µl De-Cre reaction (0.1 volumes of the chemical competent cell suspension) can be directly transformed into 200 µl chemically competent cells.*

For electrotransformation, up to 2 µl De-Cre reaction can be directly mixed with 100 µl electrocompetent cells, and transformed by using an electroporator (e.g. BIORAD *E. coli* Pulser) at 1.8-2.0 kV.

*Larger volume of De-Cre reaction must be desalted by ethanol precipitation or PCR purification column prior to electrotransformation. The desalted De-Cre reaction mix should not exceed 0.1 volumes of the electrocompetent cell suspension.*

*The cell/DNA mixture can be immediately used for electrotransformation without prior incubation on ice.*

5. Add up to 400 µl of LB media (or SOC media) per 100 µl of cell/DNA suspension immediately after the transformation (heat shock or electroporation).
6. Incubate the suspension in a 37°C shaking incubator (recovery).

*For recovery of partially deconstructed double/triple fusions, incubate the suspension in a 37°C shaking incubator for 1 to 2 hours.*

*For recovery of individual educts, incubate the suspension in a 37°C shaking incubator overnight or for at least 4 hours.*

7. Plate out the recovered cell suspension on agar containing the desired (combination of) antibiotic(s). Incubate at 37°C overnight.
8. Colonies after overnight incubation can be verified directly by restriction digestion at this stage (refer to steps 12-16).

*This is especially recommended in cases where only a single educt or partially deconstructed multifusion plasmid is desired.*

For further selection by single antibiotic challenge on a 96 well microtiter plate, continue with step 9.

*Several different single educts/partially deconstructed multifusion plasmids can be processed and selected in parallel on one 96 well microtiter plate.*

9. For 96 well analysis, inoculate four colonies each from agar plates containing a defined set of antibiotics into approx. 500 µl LB medium without antibiotics. Incubate the cell cultures in a 37°C shaking incubator for 1-2 hours.



10. While incubating the colonies, fill a 96 well microtiter plate with 150 µl antibiotic-containing LB medium or dye (positional marker) in the corresponding wells.

*Refer to Figures 12 and 13 for the arrangement of the solutions in the wells, which are used for parallel selection of single educts or partially deconstructed multifusion plasmids. The concept is that every cell suspension from a single colony needs to be challenged by all four antibiotics separately for unambiguous interpretation.*

11. Add 1 µl aliquots from the pre-incubated cell cultures (Step 9) into the corresponding wells. Incubate the 96 well microtiter plate in a 37°C shaking incubator overnight at 180-200 rpm.

*Recommended: use parafilm to wrap the plate to prevent desiccation.*

*The remainder of the pre-incubated cell cultures can be kept at 4°C in a refrigerator for further inoculations if necessary.*

12. Select transformants containing desired single educts or partially deconstructed multifusion plasmids according to the combination of dense (growth) and clear (no growth) cell cultures from each colony. Inoculate 10-20 µl from the cell cultures into 10 ml LB media with corresponding antibiotic(s). Incubate in a 37°C shaking incubator overnight.
13. The next day, centrifuge the overnight cell cultures at 4000g for 5-10 minutes. Purify plasmid from cell pellets with common plasmid miniprep kits, according to manufacturers' protocols.
14. Determine the concentrations of purified plasmid solutions by using UV absorption spectroscopy (e.g. NanoDrop™ 1000).
15. Digest 0.5-1 µg of the purified plasmid solution in a 20 µl restriction digestion (with 5-10 units endonuclease). Incubate under recommended reaction condition for approx. 2 hours.
16. Use 5-10 µl of the digestion for analytical agarose gel (0.8-1.2%) electrophoresis. Verify plasmid integrity by comparing the *de facto* restriction pattern to the *in silico* predicted restriction pattern (e.g. by using VectorNTI, Invitrogen, or any other similar program).
17. Optional: Occasionally, a deconstruction reaction is not complete but yields partially deconstructed fusions which still retain entities to be eliminated. In this case, we recommend to pick these partially deconstructed fusions containing and perform a second round of Cre deconstruction reaction (repeat steps 1-8) by using this construct as starting material.

## **D.4. Transposition protocol for pACEBac derivatives (electroporation).**

### Reagents:

Electro-chemically competent DH10VLP Factory™ cells  
(Ampicillin)  
Kanamycin  
Tetracyclin  
Gentamycin  
BluoGal  
IPTG

Approximately 5-10 ng of the sequenced pACEBac derivative or of an acceptor-donor fusion is incubated on ice (15 min) with 50-100 µl electro-competent DH10VLP Factory™ cells. Following electroporation (optimum settings need to be determined for the equipment used), cells are incubated at 37 °C for over night (8-16 hours) and plated on agar containing kanamycin (50 µg/ml), gentamycin (7 µg/ml), (ampicillin (100 µg/ml)), tetracyclin (10 µg/ml), BluoGal (100 µg/ml) and IPTG (40 µg/ml). White colonies are selected after incubation at 37 °C (18 to 24 hours). Ampicillin selection is optional. Proceed to bacmid preparation for insect cell infection (D.5.).

## **D.5. Bacmid preparation and infection of insect cells.**

Preparation of bacmid DNA, infection of insect cells and protein expression is carried out according to established protocols, e.g. O'Reilly, D.R., Miller, L.K. & Luckow, V.A. "Baculovirus expression vectors. A laboratory manual." Oxford University Press, New York - Oxford, 1994 or David W. Murhammer (ed.). Baculovirus and Insect Cell Expression Protocols, 2<sup>nd</sup> edition, Methods in Molecular Biology™ 388, Humana Press, Totowa 2007. You may also refer to Fitzgerald et al. (2006), especially pp. 1025-27 and Bieniossek et al. (2008) for protocols on insect cell culture and bacmid preparation.

Large plasmid kits (e.g. Qiagen) can be used to extract recombinant bacmids. Note that precipitation of DNA with isopropanol from cleared lysated is recommended instead of using columns.

## D.6. Determining baculoviral expression

As protein expression in the baculoviral system is driven by the late to very late polyhedrin and p10 promoters, onset of protein expression will show some lag and should be expected to start yielding significant amounts at 15 to 24 hrs post-infection. Expression will likely peak at around 40 hrs post-infection but this may vary. Therefore, rapid and reliable methods to determine peak expression are needed. Approaches in use include immunoblot analysis, protein staining techniques or metabolic labeling to chart protein accumulation. Most of these come with a delay due to the necessary handling, e.g. casting gels, preparing protein samples etc.

Expression of target genes from the VLP Factory™ bacmid can be determined by SDS-PAGE with or without subsequent Western Blot. Incorporation of tags such as 6xHis (Hexa-His) or FLAG allow purification of proteins and protein complexes. Additionally, you may equip any of your expressed proteins with a reporter protein, e.g. fluorescent proteins such as GFP or YFP. **Keep in mind that the virus itself constitutively expresses mCherry from the backbone.**

## D.7. Protocol for Influenza VLP purification

1. At the end of the expression (around 72 hours post cell division arrest, due to mCherry, the cells will have a red color similar to red wine), dispose cells in fresh Falcon tube and spin down 10 minutes @ 4000 rpm on table top centrifuge.
2. Repeat step 1.
3. Ultracentrifuge clear supernatant overnight @ 26000rpm in SW28 rotor (6 tubes of 38mL).
4. On ice, gently resuspend each pellet in 100uL PBS.
5. Prepare a 38mL ultracentrifuge tube containing 20% sucrose and around 2mL of 60% sucrose at the bottom (will form a layer called “pillow”).
6. Load 1mL of VLP pellet.
7. Run ultracentrifuge overnight @ 26000rpm in SW28 rotor (note that the tube for balance must be prepared the same way).
8. Remove 20% sucrose solution (by pipetting gently) and collect the white layer between the two sucrose solutions (interface of the pillow).

9. Load the collected layer on a discontinuous 20-60% sucrose gradient.
10. Run ultracentrifuge overnight @ 26000rpm in SW28 rotor
11. White visible band is obtain and collected.

Note 1: If removal of the sucrose is desired, the white band can be diluted in PBS, ultracentrifuged overnight @ 26000rpm in SW28 rotor, and pellet resuspended in PBS.

Note 2: All steps are performed on Ice and sucrose solutions are prepared in PBS and filtered on 0.2  $\mu\text{m}$ .

## E. Appendix

### E.1. Verifying bacmid integrity in DH10VLP Factory™ cells

Streak DH10VLP Factory™ cells (from agar stock) on LB plates containing kanamycin (50 µg/ml), tetracycline (10 µg/ml), Bluogal (100 µg/ml) or X-Gal (20 µg/ml) and IPTG (40 µg/ml) and incubate overnight at 37°C. All colonies should turn blue. Note that the dye reaction can take >24 hrs to develop completely. Working stock of bacmid-containing *E.coli* can then be prepared from individual colonies. If, after careful plating, a significant number of colonies show up white, then the stock is likely contaminated and you are advised to re-order bacmid stock (which will be provided free of charge).

### E.2. Preparing competent DH10VLP Factory™ cells

#### E.2.1 Protocol for electrocompetent cells

**Reagents required:**

Kanamycin: 50mg/ml stock solution in water (1000x)

Tetracyclin: 12 mg/ml stock solution in 70% EtOH (500x)

BluoGal: 100mg/ml stock (1000x)

IPTG: 0.5M (or 1M) stock (1000x)

Sterile 10% glycerol (mix glycerol with MilliQ water, autoclave)

Protect plates from light, store at 4°C

1. Streak out DH10VLP Factory™ cells from frozen stock or agar stab onto an agar plate with antibiotics (Kan, Tet), X-Gal or BluoGal and IPTG.

**All colonies must turn BLUE (can take 24 to 30hrs, sometimes longer if the stabs are old). If there are white colonies, the frozen stock is contaminated.**

2. Pick one to three blue colonies, prepare a 100mL preculture (make 2 glycerol stocks 30% from preculture) and inoculate 1L LB or 2xTY culture medium adding Kan and Tet as antibiotics. Shake over night at 37 °C or RT.
3. Let culture grow to OD600 = 0.5.

4. Chill 15 min on ice. Pellet by centrifuging at 4000rpm, 4°C, 15min.
5. Resuspend in 500 ml ICE COLD STERILE 10% Glycerol-solution.
6. Centrifuge at 4000rpm, 4°C, 15 min.
7. Resuspend in 250 ml ICE COLD STERILE 10% Glycerol-solution.
8. Centrifuge at 4000rpm, 4°C, 15 min.
9. Resuspend in 50 ml ICE COLD STERILE 10% Glycerol-solution.
10. Centrifuge at 4000rpm, 4°C, 15 min (in 50ml Falcon tube, table top centrifuge).
11. Resuspend in 2 ml ICE COLD STERILE 10% Glycerol-solution.
12. Prepare 100 µl aliquots (in sterile Eppendorf tubes, labeled either MBT or MBY)
13. Flashfreeze in liquid nitrogen, store at –70°C
14. Test cells by streaking out (!!) from one frozen aliquot on agar plate containing antibiotics Kan, Tet plus X-Gal or BluoGal/ IPTG.

**All colonies must be BLUE (can take 24 to 30 hrs). If there are white colonies, the competent cells are unusable and must be made anew.**

15. Test one frozen aliquot by „blank“ electroporation (1.8kV) – actual kV should be 4-4.5 or more.

You may also want to take a look at other protocols, e.g. the protocol in Sambrook and Russell's Molecular Cloning: A Laboratory Manual or the protocol for enhanced electrotransformation efficiency by Wu et al. in the Electronic Journal of Biotechnology.

## E.2.2 Protocol for chemically competent cells

Aside from the protocol provided below, there are numerous protocols for preparing competent cells for heat-shock transformation, e.g. the classical CaCl<sub>2</sub> method, the TSS protocol (originally by Chung et al., 1989, PNAS USA, Vol. 86, pp. 2172-2175), etc.

You will find various protocols online, e.g.

[http://francois.schweisguth.free.fr/protocols/TSS\\_competent\\_cell\\_preparation.pdf](http://francois.schweisguth.free.fr/protocols/TSS_competent_cell_preparation.pdf);

[http://www.personal.psu.edu/dsg11/labmanual/DNA\\_manipulations/Comp\\_bact\\_by\\_TSS.htm](http://www.personal.psu.edu/dsg11/labmanual/DNA_manipulations/Comp_bact_by_TSS.htm);

[http://openwetware.org/wiki/Preparing\\_chemically\\_competent\\_cells](http://openwetware.org/wiki/Preparing_chemically_competent_cells)

or in Sambrook and Russell.

### Preparing competent *E.coli* (TOP10) cells with the CaCl<sub>2</sub> method

1. Grow *E.coli* strain by streaking on LB agar plates; incubate o/n at 37 ° C
2. Grow a few colonies in 20ml LB at 37 ° C o/n (to OD 600=1.8)
3. Culture 13ml into 100ml freshly prepared LB (to OD 600 = 0.25)
4. Incubate at 37 ° C on a shaker (180 rpm) for 1 hour (OD600=0.5-0.6, depending on the strain)
5. Distribute the culture into 50ml centrifuge tubes and centrifuge at 8000 rpm, 4 ° C for 5 min (or, alternatively, in 50 ml Falcon tubes, spin at -4000rpm for 6 min)
6. Discard the supernatant completely; then add 10ml of ice cold 0.1M CaCl<sub>2</sub> (keep the cells on ice and resuspend gently)
7. Incubate the solution on ice for 15 min, then centrifuge at 8000 rpm at 4 ° C for 5 min (or, alternatively, in 50 ml Falcon tubes, spin at -4000rpm for 6 min)
8. Remove the supernatant completely and let the pellet air dry
9. Add 2.8ml ice cold 0.1M CaCl<sub>2</sub> and mix gently. Keep on ice till use **OR**
10. Add 1.2 ml of sterile ice cold 80% glycerol to a tube with 2.8ml cells each and mix gently; transfer 100 µl aliquots into sterile 1.5 ml Eppendorf microcentrifuge tubes (use gloves and Bunsen burner flame); keep the Eppendorfs in ice until snap-freezing the bacterial cells in liquid N<sub>2</sub>; store at -80°C. **Alternatively**, transfer 100 µl aliquots into sterile 1.5 ml Eppendorfs, add 32 µl of 80% glycerol , freeze in dry ice/EtOH bath or liquid N<sub>2</sub> and store at -80°C.

## E.3. Preparing bacterial stocks from agar stabs

We recommend preparing your personal bacterial stock from the agar stabs you received in the kit or transforming your laboratory strain of choice with the vectors (please note that for the donor vectors this needs to be a pir<sup>+</sup> strain). This is advisable since agar stabs only have a limited shelf life (3-12 months depending on storage conditions). Permanent cultures in glycerol or DMSO, on the other side, should yield viable colonies even after 1 year and beyond if properly stored.

To generate your bacterial stock for long-term storage, streak bacteria from the agar stab onto an appropriate selective plate (refer to the vector maps for acceptor and donor vectors) or plates without antibiotics (pir<sup>HC</sup> and pir<sup>LC</sup> strains; we recommend to test these strains against a panel of antibiotics to be on the safe side; no growth of colonies should be observed under conditions of antibiotic selection). Incubate the plates over night at 37°C and then proceed to prepare stocks from

individual colonies for long-term storage according to your protocol of choice (glycerol, DMSO, etc.), as described, for example, in Inoue et al. (1990), Molecular Cloning (Sambrook and Russell, 2000), Current Protocols in Molecular Biology (Ausubel et al., 1994), etc.



## E.4. VLP Factory™ vectors: maps, sequences, restriction


Note: All acceptor and donor vector sequences can be provided in electronic format. These sequences contain all relevant information such as unique restriction sites, oris, resistance markers, etc. that is also shown in the circle maps. Request your set of vector files and accompanying files from Geneva Biotech at [contact@geneva-biotech.com](mailto:contact@geneva-biotech.com).

In the next section, the acceptor and donor vectors are presented as circle maps and, in addition, the multiple cloning site (MCS) of each vector is shown featuring important unique restriction sites. Moreover, you will find, for the purposes of designing a restriction strategy, a non-exhaustive list of restriction endonucleases that cut once, twice or not at all. Additional restriction sites can be identified with any sequence analysis software, e.g. VectorNTI, ApE, etc. or by using online tools such as WebCutter 2.0 (<http://rna.lundberg.gu.se/cutter2>) or the NEB cutter V2.0 (<http://tools.neb.com/NEBcutter2/>).

Finally, for verifying vector identity, representative digestion patterns are listed in a table.

## E.4.1 Acceptor vectors

### E.4.1.1 pACEBac1: 2904 bp

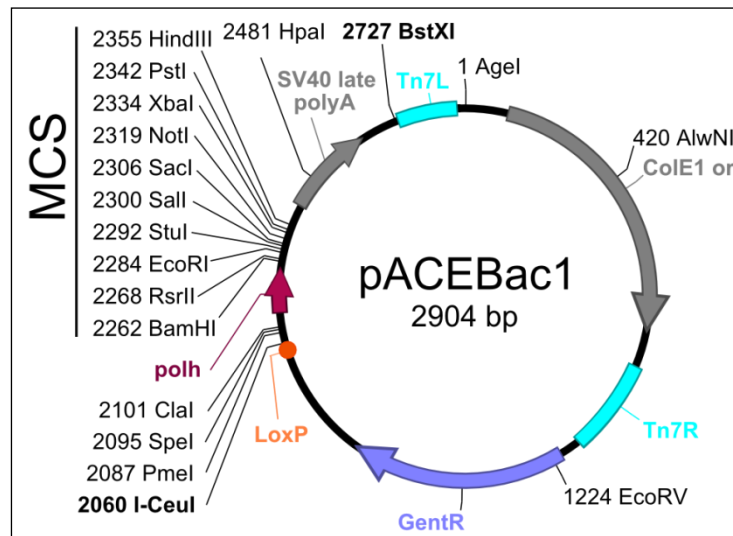

#### Multiple Cloning Site

AatI  
BamHI RsrII BssHII EcoRI StuI Sali SacI

CGGGCGCGGATCCCGGTCCGAAGCGCGCGGAATTCAAAGGCCTACGTCGACGAGCTCACTTGT

NotI BstBI XbaI PstI HindIII

CGCGGCCGCTTTTCGAATCTAGAGCCTGCAGTCTCGACAAGCTTGTCGAGAAGTACTAGAGGA

#### Enzymes that cut pACEBac1 once

2294	<b>AatI</b>	1	Age I	520	AlwNI	2127	Asel, AsnI
1629	AspI	2715	AvrII	2262	<b>BamHI</b> , BstI	2310	BanII
2619	BclI	1307	BsaAI	2102	Bsp106I, BspXI	2277	<b>BssHII</b>
2329	<b>BstBI</b>	2744	BstXI	2075	Bsu36I	2102	ClaI
2284	<b>EcoRI</b>	1226	EcoRV	890	FspI	2355	<b>HindIII</b>
2483	HpaI	2470	MfeI	1111	MscI, MluNI	2320	<b>NotI</b>
2109	NruI	2090	PmeI	2346	<b>PstI</b>	114	RcaI
2269	<b>RsrII</b>	2310	<b>SacI</b>	1169	SacII	2300	<b>Sall</b>
2370	Scal	2095	SpeI	2294	<b>StuI</b>	2334	<b>XbaI</b>

**Bold type:** restriction enzymes cutting in the MCS

#### Enzymes that cut pACEBac1 twice

2301, 2722	AccI	949, 1418	BglII	2118, 2228	BseAI, BsiMI
2470, 2569	BsmI	1169, 2320	BstZI	2118, 2228	MroI
2125, 2225	SspI				


### Enzymes that do not cut pACEBac1

Acc65I	AflII	Apal	BclI	BglI	BstEI	EcoNI	MluI
NaeI	NarI	NcoI	NdeI	NheI	NsiI	PvuI	SfiI
Smal/XmaI	SnaBI	SphI	XcmI	XhoI	XmnI		

### Sequence

5' –

accggttgacttgggtcaactgtcagaccaagtttactcatatatacttttagattgatttaaaacttcattt  
ttaattttaaaggatctaggtgaagatcctttttgataatctcatgaccaaactcccttaacgtgagttttc  
gttccactgagcgtcagacccccgtagaaaagatcaaaggatcttcttgagatccttttttctgcgcgtaat  
ctgctgcttgcaacaaaaaaaccaccgctaccagcgggtggtttgtttgccggatcaagagctaccaactct  
ttttccgaaggtaactggcttcagcagagcgcagataccaaatactgttcttctagtgtagccgtagttagg  
ccaccacttcaagaactctgtagcaccgcctacatacctcgctctgctaatacctgttaccagtggtgctgc  
cagtgggcgataagtcgtgtcttaccgggttggaactcaagacgatagttaccggataaggcgcagcggtcggg  
ctgaacggggggttcgtgcacacagcccagcttggagcgaacgacctacaccgaactgagatacctacagcg  
tgagctatgagaaagcgccacgcttcccgaagggagaaaaggcggacaggtatccggttaagcggcagggtcgg  
aacaggagagcgcacgagggagcttccagggggaaacgcctggatctttatagtcctgtcgggtttcgcca  
cctctgacttgagcgtcgatttttgtgatgctcgtcagggggcgagcctatggaaaaacgccagcaacgc  
ggcctttttacgggttcttgcccttttgcgtggccttttgcacatgttctttctgctgttatccccctgattg  
acttgggtcgctcttctgtggatgctcagatgccctgcgttaagcgggtgtgggcggacaataaagtcttaa  
actgaacaaaatagatctaaactatgacaataaagtcttaaaactagacagaatagttgtaaactgaaatcag  
tccagttatgctgtgaaaaagcatactggacttttgttatggctaaagcaaactcttcatttttctgaagtgc  
aaattgcccgtcgtatttaaagaggggctggccaagggtcatgtaaagactatattcgcggttggtgacaat  
ttaccgaacaactccgcggccgggaagccgatctcggcttgaaacgaattgttaggtggcgggtacttgggtcg  
atatcaaagtgcacatcacttcttcccgtatgcccaactttgtatagagagccactgcgggatcgtcaccgtaa  
tctgcttgacgtagatcacataagcaccaagcgcgttggcctcatgcttgaggagattgatgagcgcgggtg  
gcaatgccctgcctccggtgctcgcgggagactgcgagatcatagatatagatctcactacgcggtgctca  
aacttgggcagaacgtaagccgcgagagcgccaacaaccgcttcttggtcgaaggcagcaagcgcgatgaat  
gtcttactacggagcaagttcccagaggaatcggagtcgggtgatgttgggagtaggtggctacgtctccg  
aactcacgaccgaaaagatcaagagcagcccgcatggatttgacttggtcagggccgagcctacatgtgcga  
atgatgcccatacttgagccacctaactttgttttagggcgactgccctgctgcgtaacatcggttgctgctg  
cgtaacatcggttgctgctccataaacatcaaacatcgacccacggcgtaacgcgcttgctgcttggatgcccg  
aggcatagactgtacaaaaaaacagtcataacaagccatgaaaaccgccactgcgcccgttaccaccgctgcg  
ttcgggtcaaggttctggaccagttgctgtagcgcatacgtacttgcattacagtttacgaaccgaacaggc  
ttatgtcaactgggttcgtgccttcacccgtttccacgggtgtgcgtcacccggcaaccttgggcagcagcga  
agtcgccataacttcgtatagcatatacgaagttatctgtaactataacggtcctaaggtagcagagt  
ttaaacactagtatcgattcgcgacctaactccggaatattaatagatcatggagataattaaaaatgataacc  
atctcgcaataaataagtattttactgttttcgtaacagttttgtaataaaaaaacctataaatattccgg  
attattcataccgtcccaccatcggggcgcggatcccgggtccgaagcgcgcggaattcaaaggcctacgtcga  
cgagctcacttgctgcggccgctttcgaatctagagcctgcagtcctcgacaagcttgctcgagaagtactaga  
ggatcataatcagccataccacattttagtagaggttttacttgcgtttaaaaaacctcccacacctccccctga  
acctgaaacataaaaatgaatgcaattgttggttgaacttgtttattgcagcttataatgggttacaataaaa  
gcaatagcatcacaaatttcacaaataaagcatttttttactgcattctagttgtgggttgcctcaaaactca  
tcaatgtatcttatcatgtctggatctgatcactgcttgagcctagaagatccgggtgcttaacaaagcccgga  
aaggaagctgagttggctgctgccaccgctgagcaataactatcataaccctagggtatacccatctaatt  
ggaaccagataagtgaaatctagttccaaactattttgcatttttaattttcgtatttagcttacgacgcta  
caccaggttcccatctattttgcactcttccctaataatccttaaaaaactccatttccaccctccaggt  
tcccaactattttgcgcgccaca-3'

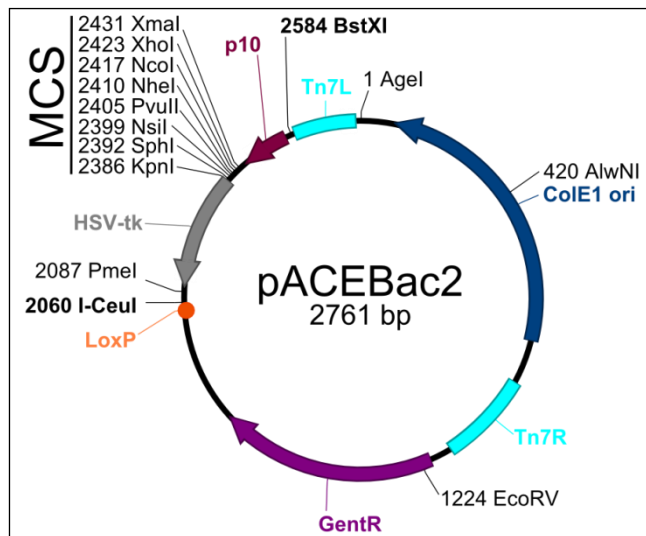

### E.4.1.2 pACEBac2: 2761 bp

#### Multiple Cloning Site

KpnI    PaeI                    NheI  
Acc65I   SphI   NsiI            PvuII

CCGTGTTTTCAGTTAGCCTCCCCATCTCCCGGTACCGCATGCTATGCATCAGCTGCTAGCA

NcoI   XhoI                    XmaI    BclI   BbsI (-)

CCATGGCTCGAGATCCCGGTGATCAAGTCTTCGTCGAGTGAT

#### Enzymes that cut pACEBac2 once

2386	<b>Acc65I</b>	1	Age I	425	AlwNI	2496	AseI, AsnI
1629	AspI	2146	BanII	2444	<b>BbsI (-)</b>	2437	<b>BclI</b>
2223	BsaI	666	BsiI	1582	BsmBI	2591	BstXI
1169	BstZI, EagI KspI	2075	Bsu36I	1226	EcoRV	2390	<b>KpnI</b>
1111	MluNI, MscI	2417	<b>NcoI</b>	2410	<b>NheI</b>	2403	<b>NsiI</b>
2090	PmeI	2407	<b>PvuII</b>	114	RcaI	1169	SacII
2433	<b>SmaI, XmaI</b>	2396	<b>SphI</b>	2423	<b>XhoI</b>		

**Bold type:** restriction enzymes cutting in the MCS    (-) cutting on the negative strand

#### Enzymes that cut pACEBac2 twice

2569, 2579	AccI	949, 1418	BglII	312, 1077	Eco57I
594, 1471	HaeII	464, 1546	HinfI		

#### Enzymes that do not cut pACEBac2

AatII	AflII	Apal	AscI	AvrII	BamHI	BclI	BglI
-------	-------	------	------	-------	-------	------	------

BssHII	Clal	Drall	EcoNI	EcoRI	HpaI	MfeI	NaeI
NarI	NdeI	NotI	NruI	PstI	PvuI	RsrII	SacI
Sall	SbfI	Scal	SfiI	SnaBI	SpeI	StuI	XbaI
XcmI	XmnI						

## Sequence

5' -

accggttgacttgggtcaactgtcagaccaagtttactcatatataacttttagattgatttaaaacttca  
 tttttaattttaaaggatctaggtgaagatcctttttgataatctcatgaccaaactcccttaacgtga  
 gttttcgttccactgagcgtcagaccccgtagaaaagatcaaaggatcttcttgagatccttttttct  
 gcgcgtaatctgctgcttgcaaaaaaaaccaccgctaccagcgggtggtttgtttgccggatcaaga  
 gctaccaactctttttccgaaggtaactggcttcagcagagcgcagataccaaatactgttcttctagt  
 gtagccgtagtttaggccaccacttcaagaactctgtagcaccgcctacatacctcgtctgctaactct  
 gttaccagtggctgctgccagtggcgataagtcgtgtcttaccgggttgactcaagacgatagttacc  
 ggataaggcgcagcggctcgggctgaacgggggggttcgtgcacacagcccagcttgagcgaacgaccta  
 caccgaactgagatacctacagcgtgagctatgagaaagcgccacgcttcccgaagggagaaaggcgga  
 caggatccggtaagcggcagggtcggaaacaggagagcgcacgaggagcgttccagggggaaacgcctg  
 gtatctttatagtcctgtcgggtttcgccacctctgacttgagcgtcgatttttgtgatgctcgtcagg  
 ggggcggagcctatggaaaaacgcagcaacgcggcctttttacgggttcctggccttttgctggccttt  
 tgctcacatgttcttttctgcttatcccttgattgacttgggtcgtcttctcgtggatgctcagatg  
 ccctgctgaagcgggtgtgggcggacaataaagtcttaaaactgaacaaaatagatctaaactatgacaa  
 taaagtcttaaaactagacagaatagttgtaaactgaaatcagtcagttatgctgtgaaaaagcatact  
 ggacttttggtatggctaaagcaaactcttcattttctgaagtgcgaattgcccgctgtattaaagagg  
 ggcgtggccaagggtcatgtaaagactatattcgcggtgtgtgacaatttaccgaacaactccgcggcc  
 gggaagccgatctcggcttgaaacgaattgttaggtggcggtacttgggtcgatatcaaagtgcatact  
 tcttcccgtatgcccactttgtatagagagccactgcgggatcgtcacccgtaactctgcttgacgtag  
 atcacataagcaccaagcgcgttggcctcatgcttgaggagattgatgagcgcgggtggcaatgcctgc  
 ctccggtgctcgcggagactgcgagatcatagatatagatctcactacgggctgctcaaacttgggc  
 agaacgtaagccgcgagagcgcgaacaacgcgttcttgggtcgaaggcagcaagcgcgatgaatgtctta  
 ctacggagcaagttcccagaggtaatcggagtcgggtgatgttgggagtaggtggctacgtctccgaac  
 tcacgaccgaaaagatcaagagcagccgcgtggttggacttgggtcagggccgagcctacatgtgcga  
 atgatgccatacttgagccacctaactttgttttagggcgactgccctgctgcgtaacatcgttgctg  
 ctgcgtaacatcgttgctgctccataacatcaaacatcgacccacggcgtaacgcgcttgctgcttgga  
 tgcccagggcatagactgtacaaaaaacagtcataacaagccatgaaaaccgcccactgcgccgttacc  
 accgctgcgttcgggtcaaggttcttgaccagttgcgtgagcgcatacgtacttgcatcagtttacg  
 aaccgaacaggccttatgtcaactgggttcgtgccttcatccgtttccacgggtgtgcgtcacccggcaac  
 cttgggcagcagcgaagtcgccataacttcgtatagcatacattatacgaagttatctgtaactataac  
 ggtcctaaggtagcaggtttaaactgacccgtagtggtatggcagggcttgccgccccgacgttggtc  
 gcgagccctgggccttcacccgaacttgggggttgggggtgggggaaaaggaagaaacgcgggctattgg  
 tcccaatgggtctcgggtgggttatcgacagagtgccagccctgggaccgaaccccgctttatgaaca  
 aacgaccaacaccgctgcgttttattctgtctttttattgccgtcatagcgcgggttccttccggtat  
 tgtctccttccgtgtttcagtttagcctccccatctcccggtaccgcatgctatgcatacagctgctagc  
 accatggctcgagatcccggtgatcaagtcttcgtcgagtgattgtaataaaaatgtaatttacagta  
 tagtatttttaattaataatacaaatgatttgataataattcttattttaactataatataattgtgttgggt  
 tgaattaaagggtccgtatactagggtatacccatctaattggaaccagataagtgaatctagttccaa  
 actattttgtcatttttaattttcgtatttagcttacgacgctacacccagttcccatctattttgtcac  
 tcttccctaataatccttaaaaactccatttccacccctccagttcccaactattttgcgcccac  
 a-3'



### Enzymes that do not cut pIDC

AatII	AflII	AlwNI	Apal	Ascl	BclI	BglI	BstZ17I
Drall	EcoNI	EcoRV	Mbol	MluI	NaeI	NarI	NdeI
NheI	NsiI	PmeI	PvuI	SbfI	SfiI	SmaI	SphI
XcmI	XhoI	XmaI	XmnI				

### Sequence

5' –

gatactagtaaaacccatgtgcctggcagataacttcgtataatgtatgctatacgaagttatgggtacc  
gcgccgcgtagaggatctgttgatcagcagttcaacctgttgatagtagtactaagctctcatgttt  
cacgtactaagctctcatgtttaacgtactaagctctcatgtttaacgaactaaacctcatggctaac  
gtactaagctctcatggctaacgtactaagctctcatgtttcacgtactaagctctcatgtttgaacaa  
taaaattaataataaatcagcaacttaaatagcctctaagggttttaagttttataagaaaaaaagaata  
tataaggcgttttaagcgttttaagggtttaacgggtgtggacaacaagccagggatgtaacgcactgaga  
agcccttagagcctctcaaagcaattttgagtgcacaggaacacttaacgggtgacagaattagcttc  
acgctgccgcaagcactcagggcgcaagggtgctaaaggaagcgggaacacgtagaaagccagtcgcga  
gaaacgggtgctgaccccgcatgaatgtcagctgggaggcagaataaatgatcatatcgtcaattattac  
ctccacggggagagcctgagcaaaactggcctcaggcatttgagaagcacacgggtcacactgcttcgggt  
agtcaataaaccggtaaacccagcaatagacataagcgggtatttaacgacctgccctgaaccgacgac  
cgggtcgaatttgctttcgaatttctgccattcatccgcttattatcacttattcaggcgtagcaacca  
ggcgtttaagggcaccaataactgccttaaaaaaattacgccccgccctgccactcatcgcagtagtgt  
tgtaattcattaagcattctgccgacatggaagccatcacaaacggcatgatgaacctgaatcgccagc  
ggcatcagcaccttgctgccttgctgataatatttgcccatgggtgaaaacggggggaagaagttgtcc  
atattggccacgttttaaatcaaaactgggtgaaactcaccagggattggctgagacgaaaaacatatcc  
tcaataaacccttttagggaaataggccagggttttcaccgtaacacgccacatcttgcaatatatgtgt  
agaaactgccggaaatcgctcggtggtattcactccagagcgatgaaaacggtttcagtttgctcatggaaa  
acgggtgtaacaagggtgaacactatcccatatcaccagctcaccgtctttcattgccatacgggaattcc  
ggatgagcattcatcaggcgggcaagaatgtgaataaaggccggataaaaacttgctgcttatttttcttt  
acgggtctttaaaaaggccgtaatatccagctgaacgggtctggttataggtacattgagcaactgactga  
aatgcctcaaaatggttctttacgatgccattgggatatatcaacgggtggtatatccagtgatttttttc  
tccatttttagcttccttagctcctgaaaatctcgataactcaaaaaatacgcccggtagtgatcttatt  
tcattatggtgaaagttggaccctcttacgtgcgatcaacgtctcattttcgccaaaagttggcccag  
atctatgtcgggtgaggagaaagaggtaatgaaatggcacctagggttatgatagttattgctcagcg  
gtggcagcagccaactcagcttccttcgggctttgttagcagccgatcttctaggctcaagcagtgga  
tcagatccagacatgataagatacattgatgagtttggaacaaaccacaactagaatgcagtgaaaaaa  
tgctttattttgtgaaattttgtgatgctattgctttattttgtaaccattataagctgcaataaacaagtt  
aacaacaacaattgcattcattttatgttttcagggttcagggggagggtgtgggagggttttttaagcaag  
taaaacctctacaaatgtggtatggctgattatgatcctctagtagtctctcgacaagcttgctcgagact  
gcaggctctagattcgaaagcggccgcgactagtgagctcgctcgacgtaggcctttgaattccgcgcgc  
ttcggaaccgggatccgcgcccgatggtgggacgggtatgaataatccggaatatttatagggtttttttat  
tacaaaactgttacgaaaacagtaaaaatacttattttatttgcgagatgggttatcattttaattatctcc  
atgatctattaataattccggagtaggtcgcaatc-3'





### Enzymes that do not cut pIDK


AatII	AflII	AgeI	Apal	ApaLI	Ascl	BamHI	BglI
Drall	EcoNI	EcoRI	EcoRV	Drall	Mbol	MluI	NdeI
PvuI	SbfI	SfiI	StuI	XbaI	XmnI		

### Sequence

5' –

gatactagtagtatacggacctttaattcaacccaacacacataatattatagttaaataagaattattatcaa  
atcatttgtatattaattaaaatactatactgtaaattacattttatttacaatcactcgcacgaagact  
tgatcacccgggatctcgcagccatggtgctagcagctgatgcatagcatgcggtaccgggagatggggg  
aggctaactgaaacacggaaggagacaataaccggaaggaacccgcgctatgacggcaataaaaagacag  
aataaaacgcacgggtgttggtcggtttgttcataaacgcggggttcggtcccagggtggtcactctgt  
cgatacccccaccgagaccccatattgggaccaataacgcccgcgtttcttccttttccccaccccaaccccc  
aagttcgggtgaaggcccgagggtcgcagccaacgctcggggcggaagccctgccatagccactacggg  
tacgttttaaaccatgtgcctggcagataacttcgtataatgtatgctatacgaagttaggtacgtac  
taagctctcatgtttcacgtactaagctctcatgtttaacgtactaagctctcatgtttaacgaactaa  
accctcatggctaacgtactaagctctcatggctaacgtactaagctctcatgtttcacgtactaagct  
ctcatgtttgaacaataaaaattaataataaatcagcaacttaaatagcctctaagggttttaagttttata  
agaaaaaaaagaatatataaggcttttaagcttttaagggttaacgggtgtggacaacaagccaggga  
tgtaacgcactgagaagcccttagagcctctcaaagcaattttcagtgacacaggaacacttaacggct  
gacagaattagcttcacgctgccgcaagcactcagggcgcaagggtgctaaaggaagcggaacacgta  
gaaagccagtcgcgcagaaacgggtgctgaccccggtatgaatgtcagctactgggctatctggacaaggga  
aaacgcaagcgcaagagaaagcaggtagcttgcagtggttacctgacgatagctagactggggcggt  
tttatggacagcaagcgaaaccggaattgccagctggggcgccctctggtaagggttggaagccctgcaa  
agtaaaactggatggctttcttgccgccaaggatctgatggcgaggggatcaagatctgatcaagagac  
aggatgaggatcggttcgcagtgattgaacaagatggattgcacgcaggttctccggccgcttgggtgga  
gaggctattcggctatgactgggcacacagacaatcggtgctctgatgccgccgtgttcgggtgctc  
agcgcagggggcgcccggttctttttgtcaagaccgacctgtccggtgccctgaatgaactgcaggacga  
ggcagcgcggctatcggtggctggccacgacgggcttccttgccgcagctgtgctcgcaggttgctactga  
agcgggaagggaactggctgctattgggcgaagtgcggggcaggatctcctgtcatctcaccttgctcc  
tgccgagaaagtatccatcatggctgatgcaatgcggcggtgcatacgttgatccggctacctgcc  
attcgaccaccaagcgaaacatcgcatcgagcgagcacgtactcggtggaagccggtccttgctgatca  
ggatgatctggacgaagagcatcaggggtcgcgccagccgaactgttcgccaggctcaaggcgcgcac  
gcccgcggcgaggatctcgtcgtgacacatggcgatgcctgcttgccgaatatcatggtggaaaatgg  
ccgcttttctggattcatcgactgtggccggctgggtgtggcgacgcgtatcaggacatagcgttggc  
taccctgatattgctgaagagcttgggcggaatgggctgaccgcttcctcgtgctttacgggtatcgc  
cgctcccgattcgcagcgcacgccttctatcgccttcttgacgagttcttctgagcgggactctgggg  
ttcgaaatgaccgaccaagcgacgccaacctgccatcacgagatttcgattccaccgccccttctat  
gaaagggttggttccggaatcggtttccgggacgcgggctggatgatcctccagcgcggggatctcatg  
ctggagttcttcgcccaccccggtatctatgtcgggtgcggagaaagaggaatgaaatggcacctagg  
tatac-3'

### E.4.2.3 pIDS: 2231 bp

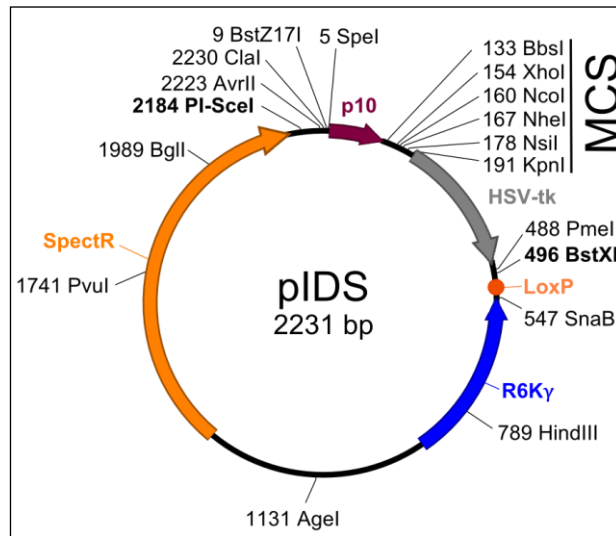

#### Multiple Cloning Site

SmaI  
BbsI      XmaI      XhoI      NcoI      NheI      NsiI  
CAATCACTCGACGAAGACTTGATCACCCGGGATCTCGAGCCATGGTGCTAGCAGCTGATGCAT

Acc65I  
SphI      KpnI  
AGCATGCGGTACC

#### Enzymes that cut pIDS once

191	<b>Acc65I, KpnI</b>	1131	AgeI	1638	ApaLI	2223	AvrII, BlnI
133	<b>BbsI</b>	1989	BglI	2087	BssHII	496	BstXI
9	BstZ17I	1081	Bsu36I	2230	ClaI	789	HindIII
2012	NaeI	160	<b>NcoI</b>	167	<b>NheI</b>	178	<b>NsiI</b>
488	PmeI	1741	PvuI	146	<b>SmaI, XmaI</b>	547	SnaBI
5	SpeI	185	<b>SphI</b>	154	<b>XhoI</b>		

**Bold type:** restriction enzymes cutting in the MCS

#### Enzymes that cut pIDS twice

82, 711	Asel, AsnI	489, 785	DraI	172, 1011	PvuII
82, 711	VspI				

#### Enzymes that do not cut pIDS

AatII	AfeI	AlwNI	Apal	Ascl	BamHI	BglII	BstBI
DraII	EcoNI	EcoRI	EcoRV	HincII	HpaI	MboI	MluI

MscI	NarI	NdeI	NotI	NruI	PstI	RsrII	SacI
Sall	SbfI	SfiI	StuI	XbaI	XcmI	XmnI	

## Sequence

5' –

cgatactagtatacggacctttaattcaacccaacacaatatattatagttaaataagaattattatca  
aatcatttgtatattaattaaaatactatactgtaaattacattttatttacaatcactcgacgaagac  
ttgatcaccggtgatctcgagccatggtgctagcagctgatgcatagcatgcggtaccgggagatgggg  
gaggctaactgaaacacggaaggagacaataaccggaaggaacccgcgctatgacggcaataaaaagaca  
gaataaaacgcacgggtgttgggtcgtttgttcataaacgcgggggttcgggtcccagggtggcactctg  
tcgataccccaccgagacccccattgggaccaataacgcccgcgtttcttcttttccccacccaacccc  
caagttcgggtgaaggcccagggtcgcagccaacgtcggggcggcaagccctgccatagccactacgg  
gtacgtttaaacccatgtgcctggcagataacttcgtataatgtatgctatacgaagtattggtacgta  
ctaagctctcatgtttcacgtactaagctctcatgtttaacgtactaagctctcatgtttaacgaacta  
aacctcatggctaacgtactaagctctcatggctaacgtactaagctctcatgtttcacgtactaagc  
tctcatgtttgaacaataaaaattaataataaatcagcaacttaaatagcctctaaggttttaagttttat  
aagaaaaaaaagaatatataaggcttttaaaagcttttaaggtttaacggttggtggacaacaagccaggg  
atgtaacgcactgagaagcccttagagcctctcaaagcaattttgagtgcacacaggaacacttaacggc  
tgacataattcagcttcacgctgccgcaagcactcagggcgcaagggtgctaaaggaagcgggaacacg  
tagaaagccagtcgcgagaaacggtgctgaccccggtgaatgtcagctgggaggcagaataaatgatc  
atatcgtcaattattacctccacggggagagcctgagcaaacctggcctcaggcatttgagaagcacacg  
gtcacactgcttcggtagtcaataaacgggtaagtagcgtatgcgctcacgcaactggtccagaacct  
tgaccgaacgcagcgggtggtaacggcgcagtgggcgttttcatggccttggtatgactgttttttgggg  
tacagtctatgcctcgggcacccaagcagcaagcgcgttacgccgtgggtcgatgtttgatgttatgga  
gcagcaacgatgttacgcagcagggcagtcgcctaaaacaaagttaaaccatcatgagggaagcgggtga  
tcgccgaagtatcgactcaactatcagaggtagttggcgtcatcgagcgccatctcgaaccgacgttgc  
tggccgtacatttgtacggctccgcagtggtggcggcctgaagccacacagtgatattgatttgctgg  
ttacggtgaccgtaaggcttgatgaaacaacgcggcgagctttgatcaacgaccttttgaaacttcgg  
cttcccctggagagagcgagattctccgcgtgtagaagtcaccattgttggtgcacgacgacatcattc  
cgtggcgttatccagctaagcgcgaactgcaatttggagaatggcagcgcaatgacattcttgcaggta  
tcttcgagccagccacgatcgacattgatctggctatcttgctgacaaaagcaagagaacatagcgttg  
ccttggtaggtccagcggcggagggaactctttgatccgggttcctgaacaggatctatttgaggcgctaa  
atgaaaccttaacgctatggaactcgccgcccgcactgggctggcgatgagcgaaatgtagtgcttacgt  
tgtcccgcatttggtacagcgcagtaaccggcaaaaatcgcgccgaaggatgtcgctgccgactgggcaa  
tggagcgcctgccggcccagtatcagccgctcatacttgaagctagacaggcttatcttggaacaagaag  
aagatcgcttggcctcgcgcgagatcagttggaagaatttgtccactacgtgaaaggcgagatcacca  
aggtagtcggcaataatgtctaacaatttcgttcaagccgacggatctatgtcgggtgcggagaaagag  
gtaatgaaatggcacctaggtat-3'

### **E.4.3. Analytical restriction digest patterns**

All VLP Factory™ acceptor and donor vectors yield characteristic restriction digest patterns when plasmid DNA is restricted with HindIII and/or HinfI or a combination of both.


<b>Vector</b>	<b>Length [bp]</b>	<b>Restriction</b>	<b>Fragment(s) [bp]</b>
<b>pACEBac1</b>	2904	HindIII	2904
		HinfI	1082, 1037, 558, 227
<b>pACEBac2</b>	2761	HindIII	2761
		HinfI	1679, 1082
<b>pIDC</b>	2312	HindIII	1697, 615
		HinfI	1125, 960, 227
		HindIII/HinfI	1100, 597, 363, 227, 25
<b>pIDK</b>	2281	HindIII	2281
		HinfI	2000, 134, 58, 52, 37
		HindIII/HinfI	1087, 913, 134, 58, 52, 37
<b>pIDS</b>	2231	HindIII	2231
		HinfI	2018, 213
		HindIII/HinfI	1413, 605, 213

#### **E.4.4 DH10VLP Factory™ baculoviral genome**

The DH10VLP Factory™ genome (size approx. 130 kb, see below schematic) is a derivative of the *Autographa californica* nucleopolyhedrovirus (AcMNPV) genome. It has been genetically engineered for improved protein production and reduced protein degradation (details above). In addition, it contains an element for accepting donor DNA into its transposition acceptor site (mini-attTn7) that concomitantly allows blue-white selection to identify successful transposition events. The bacmid possesses the F replicon from the F plasmid that keeps the plasmid copy number at 1 (single copy).

It is hosted in DH10VLP Factory™ *E.coli* cells and can be isolated from the bacteria using commercial “large construct” kits (e.g. Qiagen) or appropriate protocols for preparation of large DNA molecules (e.g. King et al., 2007). Also refer to Chapter D.5.

# VLP Factory™ Baculovirus

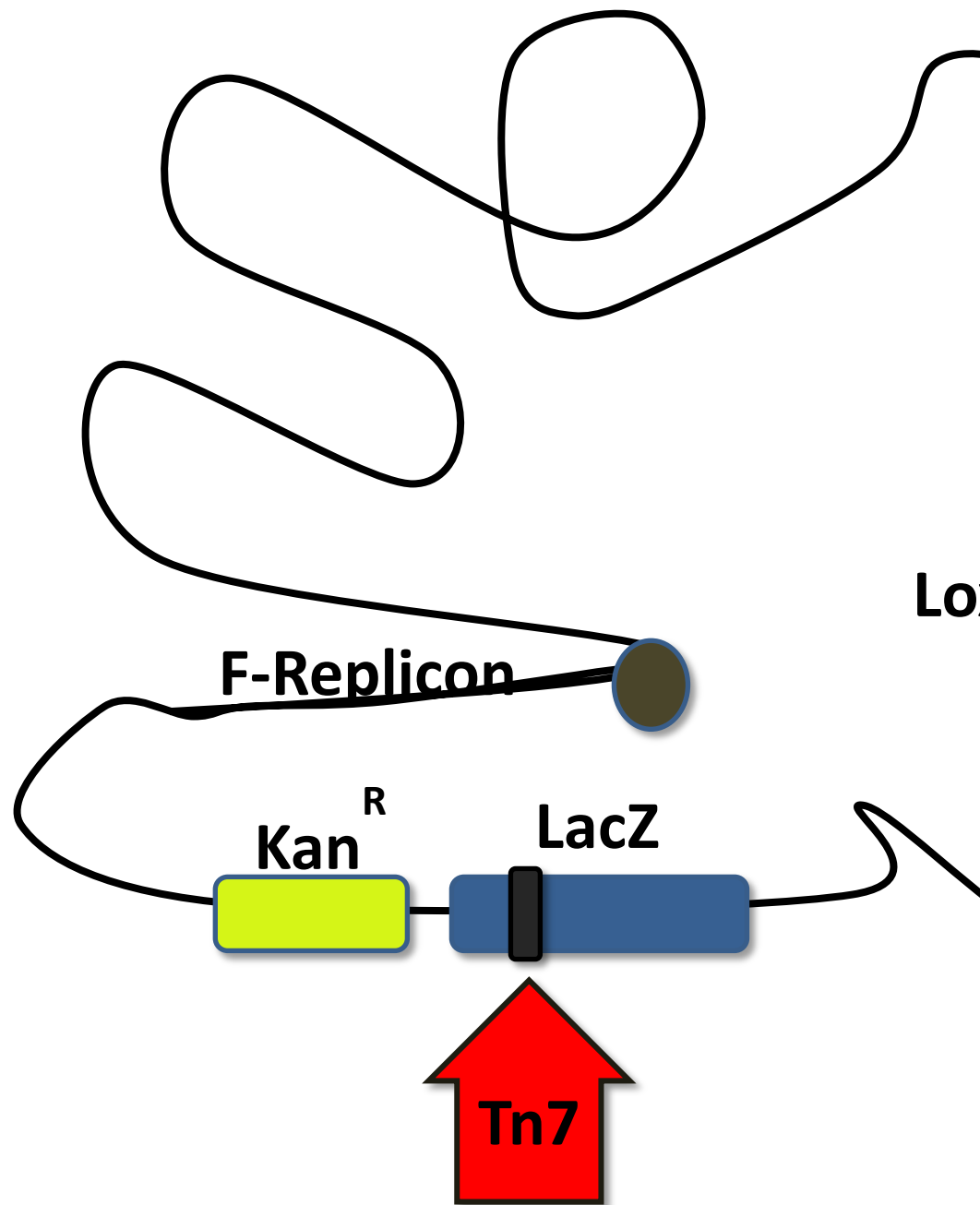

- ✓ Co-expression of Flu
- ✓ Specific Viral Gene D
- ✓ Two Separate Viral E
- ✓ Virally Encoded Fluor



## F. References

1. **Ausubel F, Brent R, Kingston RE, Moore DD, Seidman JG, Smith JA, and Struhl K** (eds., 1994). Current Protocols in Molecular Biology. John Wiley & Sons, New York, electronic version DOI 10.1002/0471142727
2. **Bieniossek C et al.** (2008). VLP Factory™ : Multigene Baculovirus-Based Eukaryotic Protein Complex Production. *Current Protocols in Protein Science* 51: 5.20.1-26
3. **Charbonnier S et al.** (2008). The social network of a cell: Recent advances in interactome mapping. *Biotechnology Annual Review* 14: 1-28.
4. **Figeys D** (2008). Mapping the human protein interactome. *Cell Research* 18: 716-724.
5. **Fitzgerald DJ et al.** (2006). Protein complex expression by using multigene baculoviral vectors. *Nature Methods* 3: 1121-1132.
6. **Hawtin RE et al.** (1995). Identification and Preliminary Characterization of a Chitinase Gene in the *Autographa californica* Nuclear Polyhedrosis Virus Genome. *Virology* 212: 673-685.
7. **Hawtin RE et al.** (1997). Liquefaction of *Autographa californica* Nucleopolyhedrovirus-Infected Insects Is Dependent on the Integrity of Virus-Encoded Chitinase and Cathepsin Genes. *Virology* 238: 243-253.
8. **Hom LG and Volkman LE** (2000). *Autographa californica* M nucleopolyhedrovirus *chiA* is required for processing of V-CATH. *Virology* 277: 178-183.
9. **Inoue H, Nojima H and Okayama H** (1990). High efficiency transformation of *Escherichia coli* with plasmids. *Gene* 96: 23-28.
10. **King L, Hitchman R, and Possee RD** (2007). Recombinant Baculovirus Isolation. In: Murhammer D (ed.): Baculovirus and Insect Cell Expression Protocols, 2<sup>nd</sup> edition, *Methods in Molecular Biology* 388: 77-93, Humana Press, Totowa, NJ.
11. **Sambrook J and Russell DW** (2000). Molecular Cloning, A Laboratory Manual, 3rd edition. Cold Spring Harbor Laboratory Press, Cold Spring Harbor, available online at <http://cshprotocols.cshlp.org>
12. **Sano K, Maeda K, Oki M, Maeda Y** (2002). Enhancement of protein expression in insect cells by a lobster tropomyosin cDNA leader sequence. *FEBS Letters* 532: 143-146
13. **Slack JM, Kuzio J and Faulkner P** (1995). Characterization of v-cath, a cathepsin L-like proteinase expressed by the baculovirus *Autographa californica* multiple nuclear polyhedrosis virus. *Journal of General Virology* 76: 1091-1098.
14. **Trowitzsch S et al.** (2010). New baculovirus expression tools for recombinant protein complex production. *Journal of Structural Biology* 172: 45-54, epub Feb 21, 2010, doi: 10.1016/j.jsb.2010.02.010
15. **Vijayachandran LS et al.** (2011). Robots, pipelines, polyproteins: Enabling multiprotein expression in prokaryotic and eukaryotic cells. *Journal of Structural Biology*
16. **Wu N et al.** (2010). Enhancing DNA electrotransformation efficiency in *Escherichia coli* DH10B electrocompetent cells. *Electronic Journal of Biotechnology* 13, doi: 10.2225/vol13-issue5-fulltext-11 (<http://ejb.ucv.cl/index.php/ejbiotechnology/article/view/v13n5-11>)



## NOTES:

## G. Purchaser Notification

### Limited Use Label License : Geneva Biotech

The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer (whether the buyer is an academic or for-profit entity).

The buyer cannot sell or otherwise transfer **(a)** this product **(b)** its components or **(c)** materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. The buyer may transfer information or materials made through the use of this product to a scientific collaborator, provided that such transfer is not for any Commercial Purpose, and that such collaborator agrees in writing **(a)** not to transfer such materials to any third party, and **(b)** to use such transferred materials and/or information solely for research and not for Commercial Purposes.

**Commercial Purposes** means any activity by a party for consideration and may include, but is not limited to: **(1)** use of the product or its components in manufacturing; **(2)** inclusion of the product as part of another product; **(3)** use of the product or its components to provide a service, information, or data; **(4)** use of the product or its components for *ex vivo* or *in vivo* therapeutic, diagnostic, prophylactic or other unauthorized commercial purposes; **(5)** use in foods, drugs, devices or cosmetics of any kind, or for consumption by or use in connection with or administration or application to humans or animals; **(6)** resale of the product or its components, whether or not such product or its components are resold for use in research, or **(7)** any use other than research use.

For products that are subject to multiple limited use label licenses, the terms of the most restrictive limited use label license shall control. Geneva Biotech will not assert a claim against the buyer of infringement of patents owned or controlled by Geneva Biotech which cover this product based upon the manufacture, use or sale of a therapeutic, clinical diagnostic, vaccine or prophylactic product developed in research by the buyer in which this product or its components was employed, provided that neither this product nor any of its components was used in the manufacture of such product.

If the purchaser does not agree with the terms or limitations of this limited use agreement, the customer can return the product at their own costs but for a full refund (excluding shipping, tax and handling fees).

For information about purchasing a license to use this product or the technology embedded in it for any use other than for research use please contact Geneva Biotech at [contact@geneva-biotech.com](mailto:contact@geneva-biotech.com)